

Contents

1. Description
 - 1.1 Background information
 - 1.2 Applications
 - 1.3 Reagent and instrument requirements
2. Protocol
3. Example of acquisition of MACSQuant Calibration Beads on a MACSQuant Analyzer

1. Description

This product is for research use only.

| | |
|-----------------------|--|
| Components | MACSQuant Calibration Beads. |
| Capacity | 3.5 mL for up to 100 tests. |
| Product format | MACSQuant Calibration Beads are supplied in aqueous solution containing 0.05% sodium azide. |
| Storage | Store bottle protected from light at +2 to +8 °C. Do not freeze. The expiration date is indicated on the vial label. |

1.1 Background information

The MACSQuant Calibration Beads are designed for routine calibration of the MACSQuant Flow Cytometers. These calibration particles are 2 and 3 µm in size and enable the adjustment of the voltage settings. In addition, the 3 µm size particles are stained with different fluorescent dyes, whose emission spectra are compatible with many commonly used fluorochromes for flow cytometry. The MACSQuant Calibration Beads are excited by light between 400 nm and 650 nm to yield emission in all fluorescent channels. This enables the successful calibration of the laser settings.

1.2 Applications

- Calibration of the scatter and fluorescence settings of the MACSQuant Flow Cytometers.

1.3 Reagent and instrument requirements

- MACSQuant Flow Cytometer, e.g., MACSQuant Analyzer 10 (# 130-096-343), MACSQuant Analyzer 16 (# 130-109-803), MACSQuant VYB (# 130-096-116), or MACSQuant X (# 130-105-100) with MACSQuantify™ Software version 2.8 or higher
- MACSQuant Running Buffer (# 130-092-747)
- Tubes, e.g., 5 mL round-bottom tubes

2. Protocol

▲ For a detailed description regarding the use of the MACSQuant Flow Cytometers refer to the respective user manual. The following guidelines are for performing an automated calibration only. Refer to the MACSQuant Flow Cytometer user manual for details concerning manual calibration.

1. Login and prime the MACSQuant Flow Cytometer. Check that the system status bar indicates 'Ready' (green).
2. Ensure that the single tube holder is correctly attached to the instrument.
3. Resuspend the MACSQuant Calibration Beads by vortexing.
4. Click on the Barcode reader icon. Position the barcode label of the MACSQuant Calibration Beads vial in front of the barcode reader.



▲ **Note:** A dialog box will appear that instructs the user to add one drop of calibration beads to a single tube.
5. Add one drop of MACSQuant Calibration Beads into an empty tube.
6. Place the tube into the single tube holder and click on run. The instrument will then proceed to automatically dilute and mix the bead suspension before performing calibration.

▲ **Note:** The system status bar will indicate 'processing sample' (blue) until completion of the task.
7. Following successful calibration the status bar will report 'MACSQuant ready; Calibration ok'; this is visually represented by a green status.

▲ **Note:** The gain and trigger settings are now preset for running experiments and can be viewed under the channel tab in the **Instrument Settings**.

3. Example of acquisition of MACSQuant Calibration Beads on a MACSQuant Analyzer


A) Forward scatter (FSC) versus side scatter (SSC)

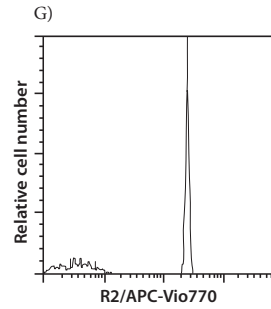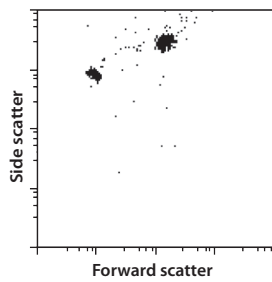

Refer to www.miltenyibiotec.com for all data sheets and protocols. Miltenyi Biotec provides technical support worldwide. Visit www.miltenyibiotec.com/local to find your nearest Miltenyi Biotec contact.

B) Fluorescence

Histograms show the fluorescence intensity of the MACSQuant Calibration Beads after excitation with the 405 nm (A), 488 nm (B, C, D, E), and 635 nm (F, G) lasers. Examples of the corresponding fluorochrome in the respective fluorescence channel are indicated in the axis legend.

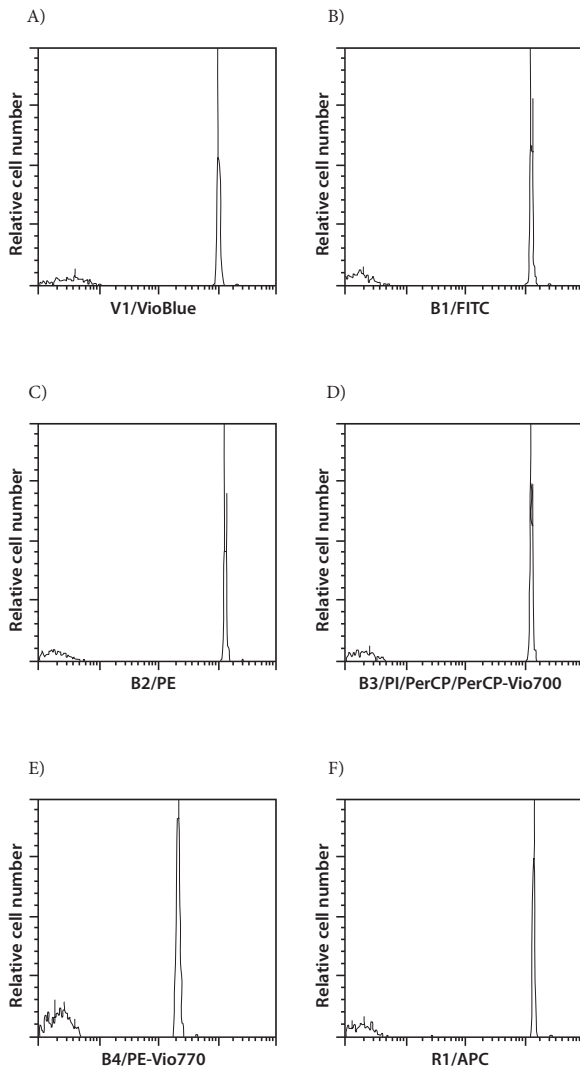

Legal notices

Limited product warranty

Miltenyi Biotec B.V. & Co. KG and/or its affiliate(s) warrant this product to be free from material defects in workmanship and materials and to conform substantially with Miltenyi Biotec's published specifications for the product at the time of order, under normal use and conditions in accordance with its applicable documentation, for a period beginning on the date of delivery of the product by Miltenyi Biotec or its authorized distributor and ending on the expiration date of the product's applicable shelf life stated on the product label, packaging or documentation (as applicable) or, in the absence thereof, ONE (1) YEAR from date of delivery ("Product Warranty"). Miltenyi Biotec's Product Warranty is provided subject to the warranty terms as set forth in Miltenyi Biotec's General Terms and Conditions for the Sale of Products and Services available on Miltenyi Biotec's website at www.miltenyibiotec.com, as in effect at the time of order ("Product Warranty"). Additional terms may apply. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING IF A PRODUCT IS SUITABLE FOR CUSTOMER'S PARTICULAR PURPOSE AND APPLICATION METHODS.

Technical information

The technical information, data, protocols, and other statements provided by Miltenyi Biotec in this document are based on information, tests, or experience which Miltenyi Biotec believes to be reliable, but the accuracy or completeness of such information is not guaranteed. Such technical information and data are intended for persons with knowledge and technical skills sufficient to assess and apply their own informed judgment to the information. Miltenyi Biotec shall not be liable for any technical or editorial errors or omissions contained herein.

All information and specifications are subject to change without prior notice. Please contact Miltenyi Biotec Technical Support or visit www.miltenyibiotec.com for the most up-to-date information on Miltenyi Biotec products.

Licenses

This product and/or its use may be covered by one or more pending or issued patents and/or may have certain limitations. Certain uses may be excluded by separate terms and conditions. Please contact your local Miltenyi Biotec representative or visit Miltenyi Biotec's website at www.miltenyibiotec.com for more information.

The purchase of this product conveys to the customer the non-transferable right to use the purchased amount of the product in research conducted by the customer (whether the customer is an academic or for-profit entity). This product may not be further sold. Additional terms and conditions (including the terms of a Limited Use Label License) may apply.

CUSTOMER'S USE OF THIS PRODUCT MAY REQUIRE ADDITIONAL LICENSES DEPENDING ON THE SPECIFIC APPLICATION. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING FOR ITSELF WHETHER IT HAS ALL APPROPRIATE LICENSES IN PLACE. Miltenyi Biotec provides no warranty that customer's use of this product does not and will not infringe intellectual property rights owned by a third party. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

Trademarks

MACSQuant, MACSQuantify, the Miltenyi Biotec logo, Vio, and VioBlue are registered trademarks or trademarks of Miltenyi Biotec and/or its affiliates in various countries worldwide.

Copyright © 2025 Miltenyi Biotec and/or its affiliates. All rights reserved.