




Miltenyi Biotec

# Flow cytometry is in their genes


REAffinity™ Recombinant Antibodies

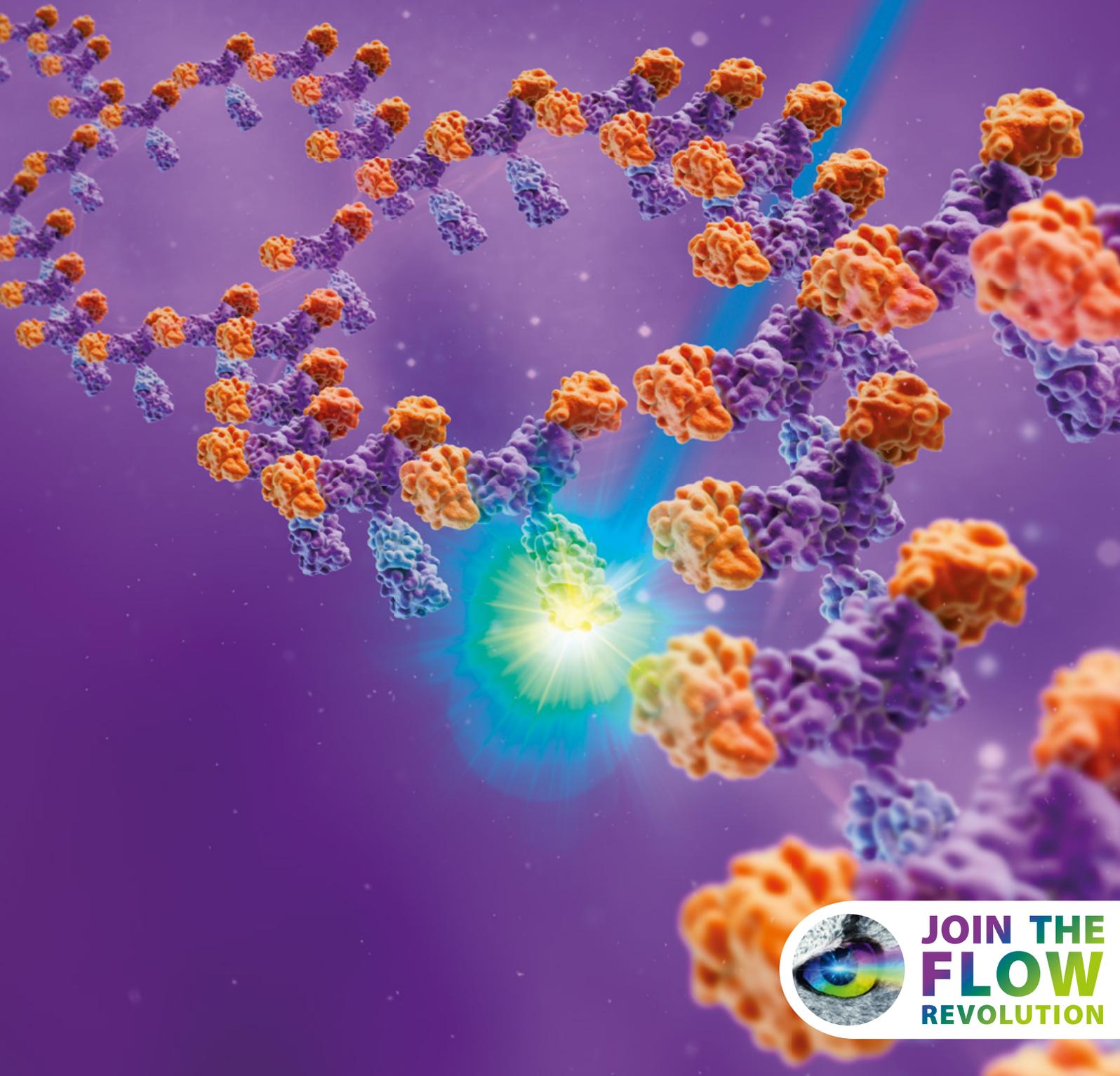

**JOIN THE  
FLOW  
REVOLUTION**

## What are REAfinity Recombinant Antibodies?

REAfinity Recombinant Antibodies are recombinantly generated antibodies that provide superior lot-to-lot consistency and purity, as compared to mouse or rat monoclonal antibodies. They have been engineered to lack any background binding. Additionally, they all have the same human IgG1 isotype, eliminating any need to include multiple isotype controls during flow analysis.

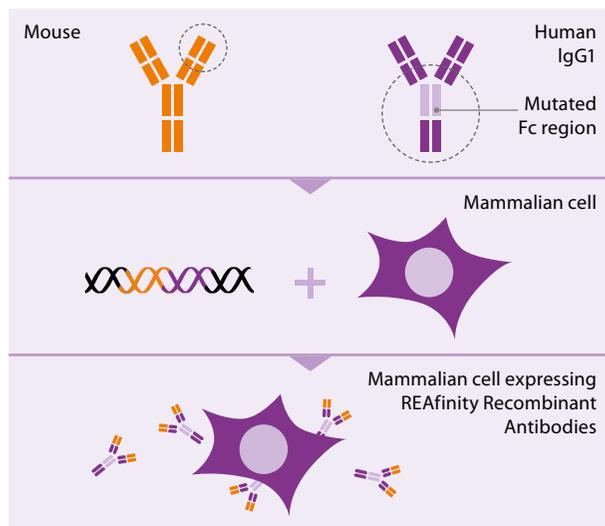

**Figure 1: Manufacturing workflow for REAfinity Recombinant Antibodies.**

REAfinity Recombinant Antibodies are generated by cloning the antigen-binding regions from traditional mouse or rat monoclonal antibodies derived from hybridomas with the human IgG1 Fc region. The two genetic regions are fused at the DNA level to create one antibody-defining genetic element (fig. 1). To eliminate background signal during analysis, the human IgG1 Fc region used for generating this fusion gene is mutated, abolishing any binding of the recombinantly engineered antibodies to FcγRs.

To guarantee maximum purity and lot-to-lot consistency, a biologically and chemically defined *in vitro* expression system is used to produce REAfinity Recombinant Antibodies. The engineered genetic sequence is expressed in a mammalian cell line that is cultured under standardized conditions.

To ensure lot-to-lot consistency, the identical cell line is used to produce all REAfinity Recombinant Antibodies, which eliminates variability due to differences in the expression system.




- ▶ High lot-to-lot consistency
- ▶ One universal isotype
- ▶ No more background signal

# Made to meet new reproducibility standards

## Get reliable results through consistent antibody quality


The manufacturing process of REAfinity Recombinant Antibodies ensures high performance consistency based on the following aspects:

- REAfinity Recombinant Antibodies are derived from a defined DNA sequence, which encodes only one type of heavy and light chain, ensuring high antibody purity (fig. 2).
- Production under biologically and chemically defined, standardized culture conditions in mammalian cells, resulting in high lot-to-lot consistency (fig. 3).
- Analytical, biochemical, and cell-based testing of every manufactured lot (fig. 3).

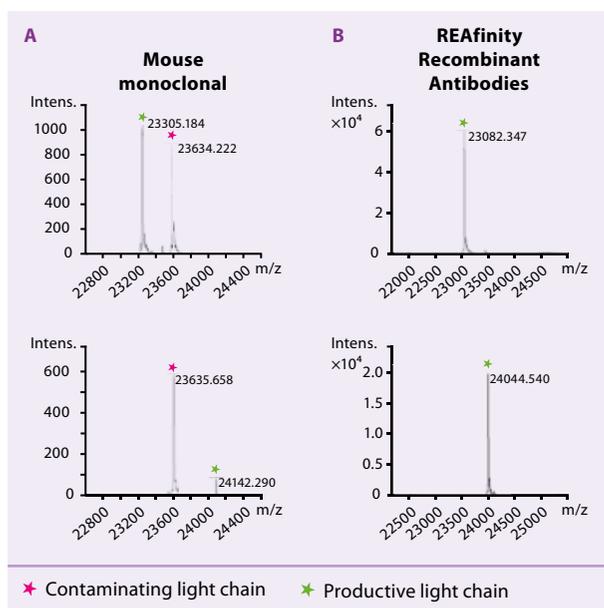

**Figure 2: Mass spectrometry analysis of purified antibodies shows that REAfinity Recombinant Antibodies are defined products, while hybridoma-generated antibodies can be a mixture.**

(A) Mass spectrometry analysis of two examples of hybridoma-generated monoclonal antibodies shows that both contain a second light chain with a molecular weight of approx. 23,635 Da. In one example bottom left, the amount of the contaminating light chain was by far exceeding the productive light chain of approx. 24,142 Da. (B) Two examples of REAfinity Recombinant Antibodies show pure light chain populations.

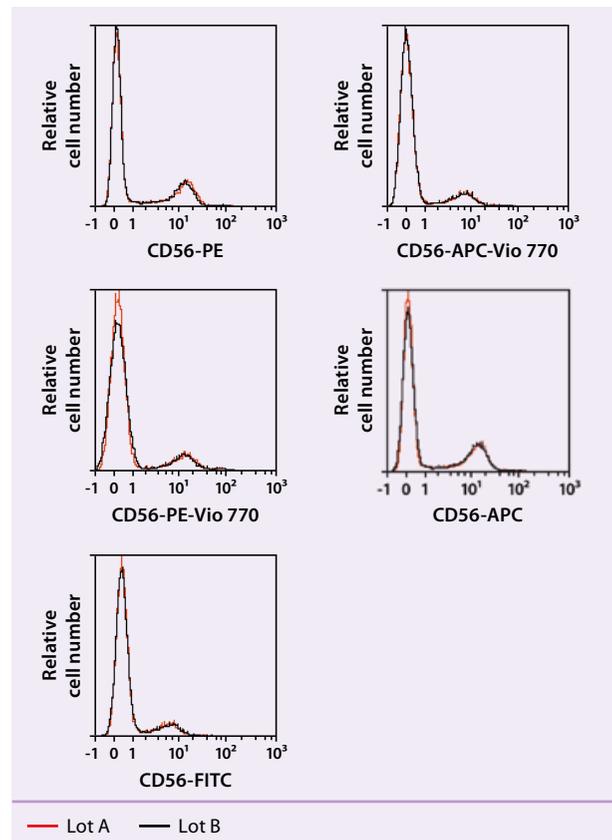


**Figure 3: Staining performance of REAfinity Recombinant Antibodies is nearly identical between different lots.** Two manufacturing lots of five different CD56 REAfinity Recombinant Antibody conjugates were compared by flow cytometry using the MACSQuant® Analyzer 10. The two histogram curves (black and red) represent the two lots. Lot-to-lot staining performance was nearly identical, when using human peripheral blood mononuclear cells (PBMCs) from a single donor.


# Hassle-free analysis – no more background signal

## Analyze what matters


All REAfinity Recombinant Antibodies contain a specifically mutated human IgG1 Fc region that abolishes their binding to Fcγ receptors (fig. 4). This allows for background-free analysis and eliminates the need for additional blocking steps such as using a FcR blocking reagent (fig. 5 and 6).

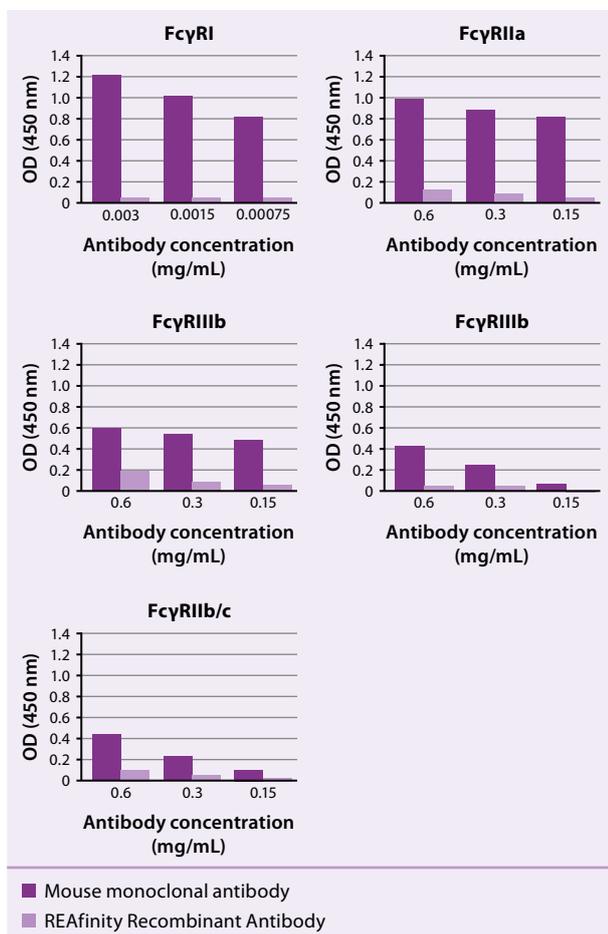

**Figure 4: Recombinantly engineered REAfinity Antibodies do not bind to Fcγ receptors.** CD144-specific hybridoma-derived mouse monoclonal antibodies (dark purple bars) bind to high affinity FcγRI (CD64) as well as low affinity FcγRIII (CD16) and FcγRII (CD32) receptors. In contrast, REAfinity Recombinant Antibodies (light purple bars) show virtually no interaction with Fcγ receptors. Binding to five cellular Fcγ receptors was compared using enzyme-linked immunosorbent assays.

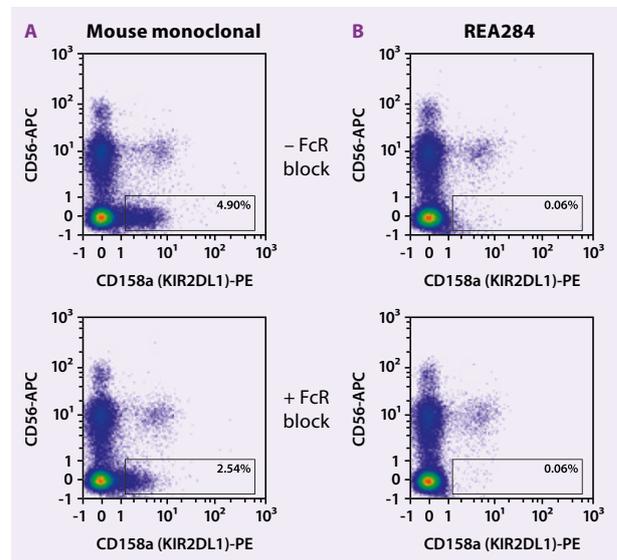

**Figure 5: Staining with REAfinity Recombinant Antibodies shows no background signal, even without FcR block.** Human PBMCs were stained with either a mouse monoclonal CD158a-PE antibody or REAfinity CD158a-PE (clone: REA284) antibody. Staining was performed without (top) and with (bottom) FcR block. The mouse monoclonal antibody binds unspecifically to CD56-negative cells (A). Staining with the REAfinity Recombinant Antibody, shows no background signal.\*

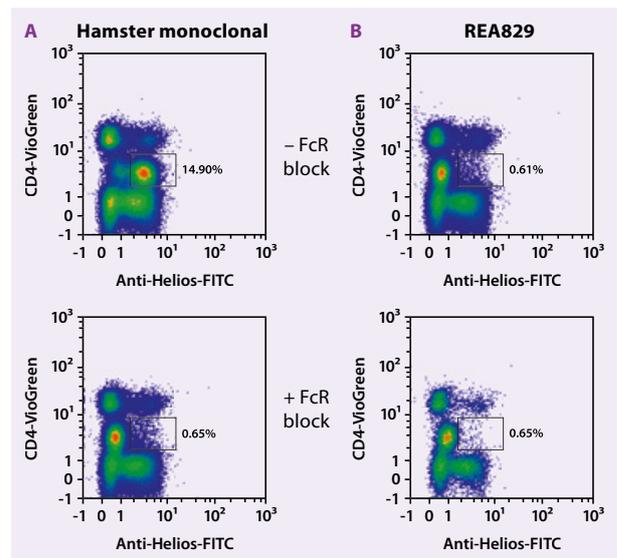


**Figure 6: Staining with REAfinity Recombinant Antibodies shows no background signal, even without FcR block.** Human PBMCs were stained with either a hamster monoclonal anti-Helios-FITC antibody or REAfinity anti-Helios-FITC (clone: REA829) antibody. Staining was performed without (top) and with (bottom) FcR block. The hamster monoclonal antibody binds unspecifically to FcγR when analyzing the CD4 intermediate monocyte population (A). Staining with the REAfinity Recombinant Antibody shows no background signal.\*

# Optimized for flow cytometry applications

## The perfect reagents for your cell analysis


REAffinity Recombinant Antibodies are designed with flow cytometry in mind. The following properties ensure their superior performance as compared to traditional hybridoma-derived monoclonal antibodies:

- Greater purity
- No background signal
- Highly optimized and standardized fluorochrome conjugation process

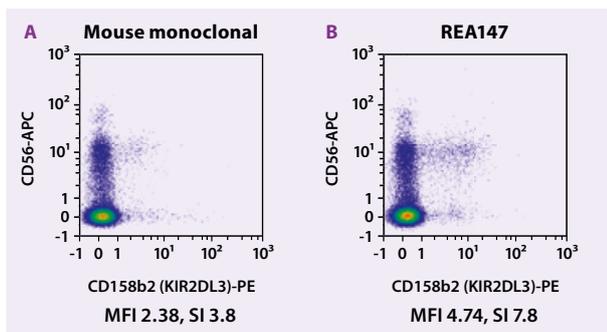


**Figure 7: Specific detection of human CD158b2<sup>+</sup> cells with clone REA147.** Human PBMCs were stained with either a PE-conjugated mouse monoclonal antibody (A) or PE-conjugated REAffinity Recombinant Antibody (B) recognizing CD158b2. Cells were also stained with CD56-APC, human (clone: REA196).\*

REAffinity Recombinant Antibody clones are available for hundreds of specificities. They are conjugated to a variety of fluorochromes to give more flexibility in setting up multiparameter cell analyses. Current options include traditional dyes, like FITC, PE, or APC, as well as Miltenyi Biotec's Vio<sup>®</sup> and Vio Bright Dyes.

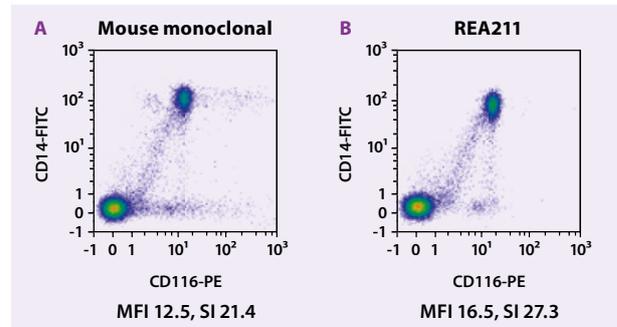

**Figure 8: Specific detection of human CD116<sup>+</sup> cells with clone REA 211.** Human PBMCs were stained with either a PE-conjugated mouse monoclonal antibody (A) or PE-conjugated REAffinity Recombinant Antibody (B) recognizing CD116. Cells were also stained with CD14-FITC.\*

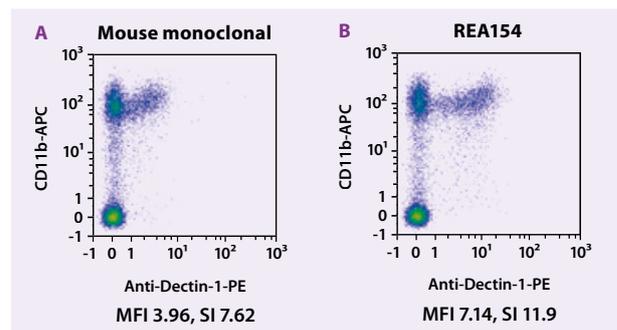

**Figure 9: Specific detection of mouse Dectin-1<sup>+</sup> cells with clone REA154.** Bone marrow cells from BALB/c mice were stained with either a PE-conjugated mouse monoclonal (A) or a PE-conjugated REAffinity Recombinant Antibody recognizing Dectin-1. Cells were also stained with CD11b-APC.\*

\* Cells were analyzed by flow cytometry on the MACSQuant Analyzer 10. Cell debris and dead cells were excluded from the analysis based on scatter signals and propidium iodide (PI) fluorescence.

# Straightforward experiment planning with only one isotype control

## Save efforts when setting up experiments

All REAfinity Recombinant Antibody clones contain the same specifically mutated human IgG1 sequence as their constant region. This offers the possibility to work with only one type of isotype control, reducing the complexity of experiment planning and saving time.

Hybridoma-derived monoclonal antibodies on the other hand, are composed of antibody isotypes derived from different species and therefore require multiple isotype controls. This complicates not only the management of reagent inventory, but also panel design when setting up experiments.

## Related products


### Isotype control antibodies

All REA clones have human IgG1 as their isotype. Therefore, only one type of isotype control, the REA Control Antibodies (clone REA293), is needed. These control antibodies are available for surface antigens (REA Control (S) Antibodies) and intracellularly expressed antigens (REA Control (I) Antibodies).

REA Control Antibodies have the same human IgG1 Fc region as the primary REAfinity Recombinant Antibodies. Thus, REA Control Antibodies are highly matched, in their structure and conjugation properties, to primary REAfinity Recombinant Antibodies. In addition, working with one isotype control reagent saves time and money.

### Compensation beads

The MACS Comp Bead Kit, anti-REA (# 130-104-693) has been developed for optimal compensation of fluorescence spillover of fluorochrome-conjugated REA clones. After staining with fluorochrome-conjugated REA clones, the MACS Comp Beads, anti-REA are used for automated or manual compensation along with the MACS Comp Beads – blank to check all the negative populations.

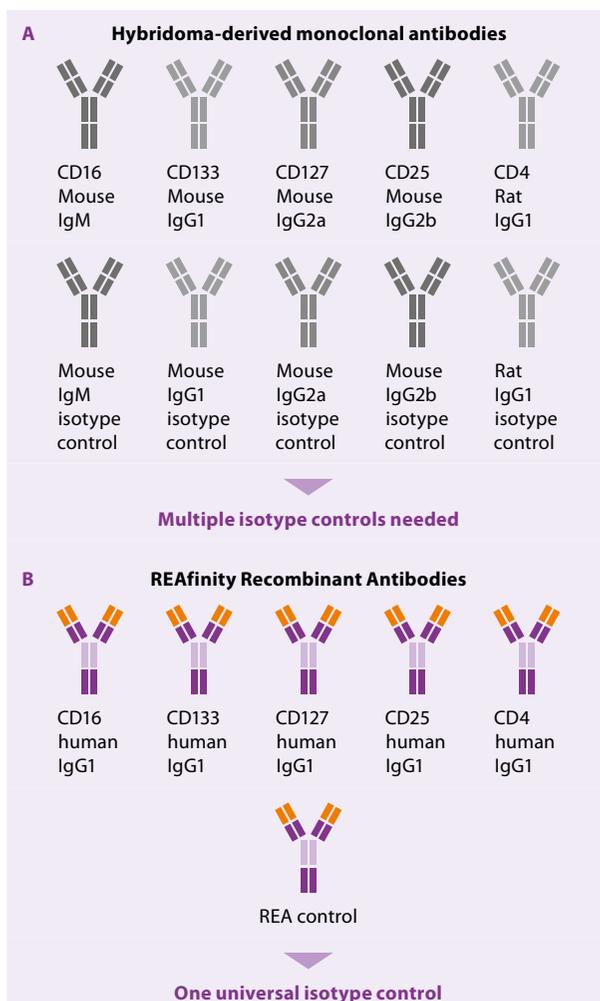

**Figure 10: REAfinity Recombinant Antibodies reduce the complexity of panel design with only one isotype control required.**



**Miltenyi Biotec**

**Germany/Austria**

Miltenyi Biotec B.V. & Co. KG  
Friedrich-Ebert-Straße 68  
51429 Bergisch Gladbach  
Germany  
Phone +49 2204 8306-0  
Fax +49 2204 85197  
[macsde@miltenyi.com](mailto:macsde@miltenyi.com)

**USA/Canada**

Miltenyi Biotec, Inc.  
2303 Lindbergh Street  
Auburn, CA 95602, USA  
Phone 800 FOR MACS  
Phone +1 866 811 4466  
Fax +1 877 591 1060  
[macsus@miltenyi.com](mailto:macsus@miltenyi.com)

**Australia**

Miltenyi Biotec  
Australia Pty. Ltd.  
Unit 11, 2 Eden Park Drive  
Macquarie Park, NSW 2113  
Australia  
Phone +61 2 8877 7400  
Fax +61 2 9889 5044  
[macsau@miltenyi.com](mailto:macsau@miltenyi.com)

**Benelux**

Miltenyi Biotec B.V.  
Sandiforddreef 17  
2333 ZZ Leiden  
The Netherlands  
[macsnl@miltenyi.com](mailto:macsnl@miltenyi.com)  
**Customer service**  
**The Netherlands**  
Phone 0800 4020120  
Fax 0800 4020100

**Customer service Belgium**

Phone 0800 94016  
Fax 0800 99626

**Customer service Luxembourg**

Phone 800 24971  
Fax 800 24984

**China**

Miltenyi Biotec Technology &  
Trading (Shanghai) Co., Ltd.  
Room A401, 4/F  
No. 1077, Zhangheng Road  
Pudong New Area  
201203 Shanghai, P.R. China  
Phone +86 21 6235 1005-0  
Fax +86 21 6235 0953  
[macs-cn@miltenyi.com.cn](mailto:macs-cn@miltenyi.com.cn)

**France**

Miltenyi Biotec SAS  
10 rue Mercœur  
75011 Paris, France  
Phone +33 1 56 98 16 16  
[macsfr@miltenyi.com](mailto:macsfr@miltenyi.com)

**Hong Kong**

Miltenyi Biotec Hong Kong Ltd.  
Unit 301, Lakeside 1  
No. 8 Science Park West Avenue  
Hong Kong Science Park  
Pak Shek Kok, New Territories  
Hong Kong  
Phone +852 3751 6698  
Fax +852 3619 5772  
[macshk@miltenyi.com.hk](mailto:macshk@miltenyi.com.hk)

**Italy**

Miltenyi Biotec S.r.l.  
Via Paolo Nanni Costa, 30  
40133 Bologna  
Italy  
Phone +39 051 6 460 411  
Fax +39 051 6 460 499  
[macsit@miltenyi.com](mailto:macsit@miltenyi.com)

**Japan**

Miltenyi Biotec K.K.  
NEX-Eitai Building 5F  
16-10 Fuyuki, Koto-ku  
Tokyo 135-0041, Japan  
Phone +81 3 5646 8910  
Fax +81 3 5646 8911  
[macsjp@miltenyi.com](mailto:macsjp@miltenyi.com)

**Nordics and Baltics**

Miltenyi Biotec Norden AB  
Medicon Village  
Scheeletorget 1  
223 81 Lund  
Sweden  
[macsse@miltenyi.com](mailto:macsse@miltenyi.com)  
**Customer service Sweden**  
Phone 0200 111 800  
Fax +46 280 72 99

**Customer service Denmark**

Phone 80 20 30 10  
Fax +46 46 280 72 99

**Customer service**

**Norway, Finland, Iceland,  
and Baltic countries**  
Phone +46 46 280 72 80  
Fax +46 46 280 72 99

**Singapore**

Miltenyi Biotec Asia Pacific Pte Ltd.  
438B Alexandra Road, Block B  
Alexandra Technopark  
#06-01  
Singapore 119968  
Phone +65 6238 8183  
Fax +65 6238 0302  
[maccs-g@miltenyi.com](mailto:maccs-g@miltenyi.com)

**South Korea**

Miltenyi Biotec Korea Co., Ltd.  
Arigi Bldg. 8F  
562 Nonhyeon-ro  
Gangnam-gu  
Seoul 06136, South Korea  
Phone +82 2 555 1988  
Fax +82 2 555 8890  
[macskr@miltenyi.com](mailto:macskr@miltenyi.com)

**Spain**

Miltenyi Biotec S.L.  
C/Luis Buñuel 2  
Ciudad de la Imagen  
28223 Pozuelo de Alarcón (Madrid)  
Spain  
Phone +34 91 512 12 90  
Fax +34 91 512 12 91  
[macs-es@miltenyi.com](mailto:macs-es@miltenyi.com)

**Switzerland**

Miltenyi Biotec Swiss AG  
Gibelinstrasse 27  
4500 Solothurn  
Switzerland  
Phone +41 32 623 08 47  
Fax +49 2204 85197  
[macsch@miltenyi.com](mailto:macsch@miltenyi.com)

**United Kingdom**

Miltenyi Biotec Ltd.  
Almac House, Church Lane  
Bisley, Surrey GU24 9DR, UK  
Phone +44 1483 799 800  
Fax +44 1483 799 811  
[macsuk@miltenyi.com](mailto:macsuk@miltenyi.com)

[www.miltenyibiotec.com](http://www.miltenyibiotec.com)

Miltenyi Biotec provides products and services worldwide. Visit [www.miltenyibiotec.com/local](http://www.miltenyibiotec.com/local) to find your nearest Miltenyi Biotec contact.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research use only and not for therapeutic or diagnostic use. MACSQuant, the Miltenyi Biotec logo and REAfinity are registered trademarks or trademarks of Miltenyi Biotec and/or its affiliates in various countries worldwide. Copyright © 2022 Miltenyi Biotec and/or its affiliates. All rights reserved.