

## Contents

1. Description
  - 1.1 Principle of the Treg Expansion Kit
  - 1.2 Background information
  - 1.3 Applications
  - 1.4 Reagent and instrument requirements
2. Protocol
  - 2.1 Sample preparation
  - 2.2 Preparation of CD3/CD28 MACSiBead™ Particles
  - 2.3 Preparation of cells and expansion
  - 2.4 Restimulation of Treg cells
  - 2.5 Removal of MACSiBead™ Particles
3. Examples of Treg cell expansion using the Treg Expansion Kit

## 1. Description

<b>Components</b>	2 mL CD3/CD28 MACSiBead™ Particles, cell culture grade, corresponding to $1.2 \times 10^8$ MACSiBead Particles pre-loaded with CD3 and CD28 antibodies.
<b>Product format</b>	All components are supplied in azide-free buffer; MACSiBead Particles contain stabilizer.
<b>Storage</b>	Store protected from light at +2 to +8 °C. Do not freeze. The expiration date is indicated on the vial label.

### 1.1 Principle of the Treg Expansion Kit

The Treg Expansion Kit, mouse is based on MACSiBead Particles pre-loaded with CD3 and CD28 antibodies. Best expansion of Treg cells is accomplished by using pre-loaded MACSiBead Particles and Treg cells at a bead-to-cell ratio of 3:1 and recombinant interleukin 2 (rIL-2) with a concentration of 2000 U/mL.

### 1.2 Background information

Regulatory T cells (Treg cells) have been described to be hypoproliferative in response to polyclonal stimulation and interleukin 2 (IL-2) *in vitro*. Thus, *in vitro* expansion of Treg cells often results in low expansion rates or low frequencies of FoxP3<sup>+</sup> Treg cells due to overgrowth by conventional T cells or loss of FoxP3 expression.

The kit is designed to efficiently expand Treg cells and to maintain FoxP3 expression after isolation with the CD4<sup>+</sup>CD25<sup>+</sup> Regulatory T Cell Isolation Kit from single-cell suspensions of mouse spleen.

Expanded Treg cells can be used for any downstream application, such as adoptive Treg cell transfer, cytokine analysis, gene expression analysis, and suppression assays.

MACSiBead Particles show no autofluorescence and normally do not need to be removed prior to flow cytometric analysis. However, if desired, removal of MACSiBead Particles is easily achieved by using the MACSiMAG™ Separator (refer to 2.5)

### 1.3 Applications

- Expansion of mouse Treg cells after isolation with the CD4<sup>+</sup>CD25<sup>+</sup> Regulatory T Cell Isolation Kit, mouse.

### 1.4 Reagent and instrument requirements

- Medium: RPMI 1640 with stable glutamine supplemented with 10% fetal bovine serum (FBS), 1% sodium pyruvate, 10 mM HEPES, and 2000 U/mL rIL-2.
  - ▲ **Note:** 2-Mercaptoethanol (0.01 mM) can be added to preserve cell viability in case of rapid cell growth.
- (Optional) CD4<sup>+</sup>CD25<sup>+</sup> Regulatory T Cell Isolation Kit, mouse (# 130-091-041).
- Recombinant interleukin 2 (rIL-2), e.g., Mouse IL-2 IS, research grade (10 µg, # 130-120-330).
- Humidified incubator.
- 96-well flat bottom plates.
- (Optional) MACSiMAG Separator (# 130-092-168) for removal of MACSiBead Particles after Treg cell expansion prior to downstream experiments.
  - ▲ **Note:** Do not remove MACSiBead Particles by using MACS® Columns and MiniMACS™, MidiMACS™, VarioMACS™, SuperMACS™, autoMACS® or autoMACS Pro Separators.
- (Optional) Fluorochrome-conjugated antibodies for flow cytometric analysis of Treg cells, for example, the Treg Analysis Kit, with FoxP3-Vio® R667, anti-mouse, REAfinity™ (# 130-120-671) or the Treg Analysis Kit, with FoxP3-PE, anti-mouse, REAfinity (# 130-120-674), including CD4 and CD25 antibodies, FoxP3 antibodies, FoxP3 Staining Buffer Set, and an optimized protocol. For more information about antibodies refer to [www.miltenyibiotec.com/antibodies](http://www.miltenyibiotec.com/antibodies).

## 2. Protocol

▲ All steps in the protocol have to be performed under sterile conditions.

▲ The protocol is intended for the stimulation of  $1 \times 10^6$  Treg cells.

### 2.1 Sample preparation

When working with lymphoid organs, non-lymphoid tissues, or peripheral blood, prepare a single-cell suspension using manual methods or the gentleMACS™ Dissociator.

For details refer to the protocols section at [www.miltenyibiotec.com/protocols](http://www.miltenyibiotec.com/protocols).

### 2.2 Preparation of CD3/CD28 MACSiBead™ Particles

1. Resuspend CD3/CD28 MACSiBead™ Particles thoroughly and transfer 50  $\mu$ L to a suitable tube.

▲ **Note:** The concentration of the Treg Expansion Kit is  $6 \times 10^7$  MACSiBead Particles per mL.

2. Add 300–600  $\mu$ L of culture medium and centrifuge at  $300 \times g$  for 5 minutes. Aspirate supernatant completely.

3. Resuspend CD3/CD28 MACSiBead Particles in 500  $\mu$ L medium (concentration of  $6 \times 10^6$  beads/mL). The reagent is ready to use.

### 2.3 Preparation of cells and expansion

▲ Start with Treg cells isolated with the CD4<sup>+</sup>CD25<sup>+</sup> Regulatory T Cell Isolation Kit, mouse. For details concerning Treg isolation refer to the respective data sheet.

1. Determine the concentration and the total number of Treg cells.  $1 \times 10^5$  cells per well are needed.

2. Transfer required volumes of cell suspension to suitable tubes.

3. Add 5–10 volumes culture medium to the cells and centrifuge at  $300 \times g$  for 10 minutes. Aspirate supernatant completely.

4. Resuspend cells at a concentration of  $2 \times 10^6$  cells/mL of medium including 2000 U/mL rIL-2. Pipette 50  $\mu$ L in a well of a 96-well flat bottom plate (day 0).

5. Add 50  $\mu$ L of CD3/CD28 MACSiBead Particles in every well.

6. At day 1, add 100  $\mu$ L media including 2000 U/mL rIL-2.

7. At day 3–5, according to media usage either split cells or aspirate 100  $\mu$ L medium and add 100  $\mu$ L medium including 2000 U/mL rIL-2.

8. For restimulation proceed to 2.4.

### 2.4 Restimulation of Treg cells

▲ For restimulation of Treg cells after 6–7 days of culture, CD3/CD28 MACSiBead Particles may be removed. Please refer to 2.5 Removal of MACSiBead Particles.

▲ **Note:** Restimulation is performed with a bead-to-cell ratio of 1:1.

Please follow steps 1–8 in section 2.3 Preparation of cells and expansion. Restimulation may also be performed in 96-well plates with a ratio of one CD3/CD28 MACSiBead Particle per Treg cell.

### 2.5 Removal of MACSiBead™ Particles

▲ Removal of MACSiBead™ Particles may be required before magnetic separation of cells with MACS MicroBeads or before restimulation with different agents or antigens.

1. Harvest cells and transfer to a 5 mL, 15 mL, or 50 mL tube and wash once with buffer.

2. Resuspend cells in buffer at a density of up to  $2 \times 10^7$  cells per 1 mL and vortex thoroughly.

3. Place the tube in the magnetic field of the MACSiMAG Separator.

▲ **Note:** Use tube rack to insert 5 mL tube into the magnetic field of the separator. For details see MACSiMAG Separator data sheet.

4. Allow the MACSiBead Particles to adhere to the wall of the tube:

5 mL tubes:	2 minutes
15 mL or 50 mL tubes:	4 minutes

5. Retaining the tube in the magnet, carefully remove the supernatant containing the MACSiBead-depleted cells and place in a new tube.

6. Remove the tube from the separator and add buffer to the same volume as before.

7. Vortex sample, replace tube in the MACSiMAG Separator and repeat steps 4–5.


8. Collected cells can now be further processed as required.

### 3. Examples of Treg cell expansion using the Treg Expansion Kit


A single-cell suspension from mouse spleen was prepared using the program `m_spleen_01` on the gentleMACS Dissociator.

Mouse Treg cells were isolated from this single-cell suspension using the CD4<sup>+</sup>CD25<sup>+</sup> Regulatory T Cell Isolation Kit and expanded for 13 days with the Treg Expansion Kit according to the protocol. Treg cells before (A) and after expansion (B) were fluorescently stained with the Treg Detection Kit (CD4/CD25/FoxP3) (APC) (# 130-094-164) and analyzed by flow cytometry using the MACSQuant<sup>®</sup> Analyzer. Expansion of Treg cells (C) was calculated from FoxP3-expressing cells on day 0, 7, and day 13, respectively.


#### A) Treg cells before expansion

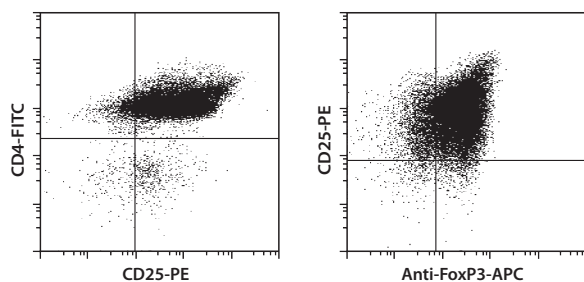

#### B) Expanded Treg cells

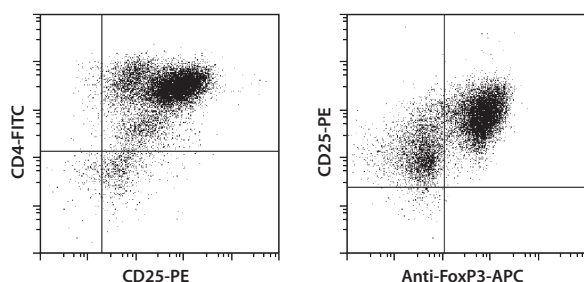

#### C) Expansion of Treg cells

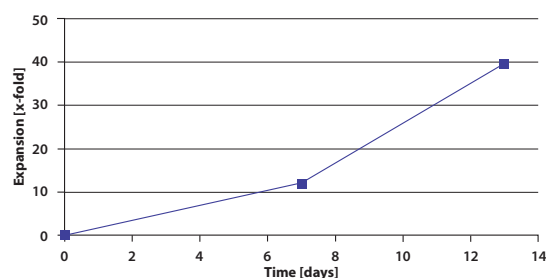

Refer to [www.miltenyibiotec.com](http://www.miltenyibiotec.com) for all data sheets and protocols. Miltenyi Biotec provides technical support worldwide. Visit [www.miltenyibiotec.com](http://www.miltenyibiotec.com) for local Miltenyi Biotec Technical Support contact information.

### Legal notices

#### Limited product warranty

Miltenyi Biotec B.V. & Co. KG and/or its affiliate(s) warrant this product to be free from material defects in workmanship and materials and to conform substantially with Miltenyi Biotec's published specifications for the product at the time of order, under normal use and conditions in accordance with its applicable documentation, for a period beginning on the date of delivery of the product by Miltenyi Biotec or its authorized distributor and ending on the expiration date of the product's applicable shelf life stated on the product label, packaging or documentation (as applicable) or, in the absence thereof, ONE (1) YEAR from date of delivery ("Product Warranty"). Miltenyi Biotec's Product Warranty is provided subject to the warranty terms as set forth in Miltenyi Biotec's General Terms and Conditions for the Sale of Products and Services available on Miltenyi Biotec's website at [www.miltenyibiotec.com](http://www.miltenyibiotec.com), as in effect at the time of order ("Product Warranty"). Additional terms may apply. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING IF A PRODUCT IS SUITABLE FOR CUSTOMER'S PARTICULAR PURPOSE AND APPLICATION METHODS.

#### Technical information

The technical information, data, protocols, and other statements provided by Miltenyi Biotec in this document are based on information, tests, or experience which Miltenyi Biotec believes to be reliable, but the accuracy or completeness of such information is not guaranteed. Such technical information and data are intended for persons with knowledge and technical skills sufficient to assess and apply their own informed judgment to the information. Miltenyi Biotec shall not be liable for any technical or editorial errors or omissions contained herein.

All information and specifications are subject to change without prior notice. Please contact Miltenyi Biotec Technical Support or visit [www.miltenyibiotec.com](http://www.miltenyibiotec.com) for the most up-to-date information on Miltenyi Biotec products.

#### Licenses

This product and/or its use may be covered by one or more pending or issued patents and/or may have certain limitations. Certain uses may be excluded by separate terms and conditions. Please contact your local Miltenyi Biotec representative or visit Miltenyi Biotec's website at [www.miltenyibiotec.com](http://www.miltenyibiotec.com) for more information.

The purchase of this product conveys to the customer the non-transferable right to use the purchased amount of the product in research conducted by the customer (whether the customer is an academic or for-profit entity). This product may not be further sold. Additional terms and conditions (including the terms of a Limited Use Label License) may apply.

CUSTOMER'S USE OF THIS PRODUCT MAY REQUIRE ADDITIONAL LICENSES DEPENDING ON THE SPECIFIC APPLICATION. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING FOR ITSELF WHETHER IT HAS ALL APPROPRIATE LICENSES IN PLACE. Miltenyi Biotec provides no warranty that customer's use of this product does not and will not infringe intellectual property rights owned by a third party. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

#### Trademarks

autoMACS, gentleMACS, MACS, MACSiBead, MACSiMAG, MACSQuant, MidiMACS, the Miltenyi Biotec logo, MiniMACS, SuperMACS, REAfinity, Vio, and VarioMACS are registered trademarks or trademarks of Miltenyi Biotec and/or its affiliates in various countries worldwide. All other trademarks mentioned in this publication are the property of their respective owners and are used for identification purposes only.

Copyright © 2026 Miltenyi Biotec and/or its affiliates. All rights reserved.