

Index

1. Description

- 1.1 Principle of MACS[®] separation
- 1.2 Background and product applications
- 1.3 Reagent and instrument requirements

2. Protocol

- 2.1 Sample preparation
- 2.2 Magnetic labeling of non-T cells
- 2.3 Magnetic separation:
Depletion of non-T cells
- 2.4 Magnetic labeling of TCR γ/δ^+ T cells
- 2.5 Magnetic separation:
Positive selection of TCR γ/δ^+ T cells

3. Example of a separation using the TCR γ/δ^+ T Cell Isolation Kit

4. Reference

1. Description

Components	1 mL Non-T Cell Depletion Cocktail, mouse: Cocktail of MicroBeads conjugated to monoclonal antibodies against mouse CD45R (B220; isotype: rat IgG2a), and CD11b (Mac-1 α-chain; isotype: rat IgG2b) and a biotin-conjugated monoclonal antibody against mouse TCRγ/δ (isotype: hamster IgG2). 1 mL Anti-Biotin MicroBeads: MicroBeads conjugated to monoclonal anti-biotin antibody (isotype: mouse IgG1).
Size	For 2×10^9 nucleated cells.
Product format	The Depletion Cocktail is supplied in a solution containing stabilizer and 0.05% sodium azide. The Anti-Biotin MicroBeads are supplied as a suspension containing 0.05% sodium azide.
Storage	Store protected from light at 4–8 °C. Do not freeze. The expiration date is indicated on the vial label.

1.1 Principle of MACS[®] separation


The isolation of mouse TCR γ/δ^+ T cells is performed in a two-step procedure. To pre-enrich the target cells, non-T cells are magnetically labeled with a cocktail of CD45R(B220) and CD11b antibodies conjugated to MicroBeads. Concomitantly, TCR γ/δ^+ T cells are labeled with Anti-TCR γ/δ -Biotin included in the cocktail. The magnetically labeled, unwanted cells are subsequently depleted by separation over a MACS[®] Column.

In the second step, TCR γ/δ^+ T cells are indirectly magnetically labeled with Anti-Biotin MicroBeads and isolated by positive selection from the pre-enriched T cell fraction. The magnetically labeled TCR γ/δ^+ T cells

are retained on the column and eluted after removal of the column from the magnetic field.

1.2 Background and product applications

The TCR γ/δ^+ T Cell Isolation Kit has been developed for the isolation of TCR γ/δ^+ T cells from single-cell suspensions of spleen, lymph nodes, or thymus. T cells bearing γ/δ TCRs are a phenotypically and functionally conserved lineage of T cells, which often constitutes the major T cell population in epithelial-rich tissues, e.g. intestine, skin, lung, and reproductive tract. In blood and lymphoid tissues, they are found at low frequency. Depending on their tissue distribution and local microenvironment, TCR γ/δ^+ T cells are involved in immune surveillance against tumors, wound healing and pro- and anti-inflammatory responses against pathogens.¹

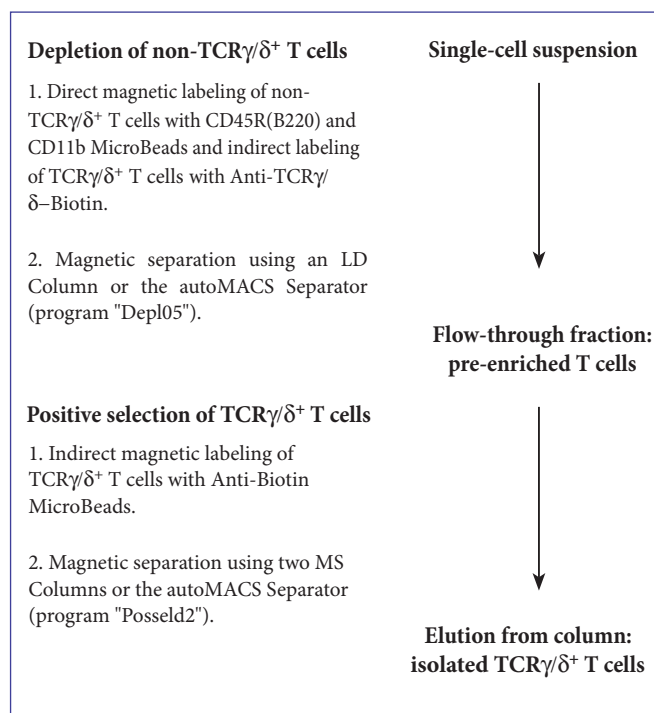

Examples of applications

Isolation of TCR γ/δ^+ T cells from single-cell suspensions of lymphoid organs for:

- phenotypical and functional analysis;
- studies on cytokine expression and receptor signaling;
- adoptive transfer or co-culture experiments.

1.3 Reagent and instrument requirements

- Buffer (degassed): Prepare a solution containing PBS (phosphate buffered saline) pH 7.2, 0.5% BSA and 2 mM EDTA by diluting MACS BSA Stock Solution (# 130-091-376) 1:20 with autoMACS[™] Rinsing Solution (# 130-091-222). Keep buffer cold (4–8 °C).

▲ **Note:** EDTA can be replaced by other supplements such as anticoagulant citrate dextrose formula-A (ACD-A) or citrate phosphate dextrose (CPD). BSA can be replaced by other proteins such as gelatine, mouse serum, or fetal calf serum. Buffers or media containing Ca^{2+} or Mg^{2+} are not recommended for use.

- MACS Columns and MACS Separators: Depletion of non-TCR γ/δ^+ T cells is performed on an LD Column. The subsequent positive selection of TCR γ/δ^+ T cells is performed on two MS Columns. Depletion and positive selection can also be performed by using the autoMACS Separator.

Column	max. number of labeled cells	max. number of total cells	Separator
Depletion			
LD	10^8	5×10^8	MidiMACS, QuadroMACS, VarioMACS, SuperMACS
Positive selection			
MS	10^7	2×10^8	MiniMACS, OctoMACS, VarioMACS, SuperMACS
Positive selection or depletion			
autoMACS	2×10^8	4×10^9	autoMACS

▲ **Note:** Column adapters are required to insert certain columns into VarioMACS™ Separator or SuperMACS™ Separator. For details, see MACS Separator data sheets.

- (Optional) Fluorochrome-conjugated antibodies for flow cytometric analysis, e.g. CD3 and Anti-TCR γ/δ .
- (Optional) Propidium iodide (PI) or 7-AAD for flow cytometric exclusion of dead cells.
- (Optional) Pre-Separation Filters (# 130-041-407) to remove cell clumps.

2. Protocol

2.1 Sample preparation

Prepare a single-cell suspension from lymphoid organs using standard methods.

▲ **Note:** Dead cells may bind non-specifically to MACS™ MicroBeads. In case of high numbers of dead cells, removal of dead cells by density gradient centrifugation or the Dead Cell Removal Kit (# 130-090-101) is recommended.



2.2 Magnetic labeling of non-T cells

- Work fast, keep cells cold and use pre-cooled solutions. This will prevent capping of antibodies on the cell surface and non-specific cell labeling.
- Volumes for magnetic labeling given below are for up to 10^8 leukocytes. When working with fewer than 10^8 cells, use the same volumes as indicated. When working with higher cell numbers, scale up all reagent volumes and total volumes accordingly (e.g. for 2×10^8 leukocytes use twice the volume of all indicated reagent volumes and total volumes).
- For optimal performance it is important to obtain a single-cell suspension before magnetic separation. Pass cells through 30 μm nylon mesh (Pre-Separation Filters # 130-041-407) to remove cell clumps which may clog the column.

1. Determine the number of leukocytes.
2. Centrifuge cells at $300 \times g$ for 10 minutes. Pipette off supernatant completely.
3. Resuspend cell pellet in 450 μL of buffer per 10^8 total cells.
4. Add 50 μL of **Non-T Cell Depletion Cocktail, mouse** per 10^8 total cells.
5. Mix well and incubate for 15 minutes at $4-8^\circ\text{C}$.
6. Wash cells by adding $10-20 \times$ labeling volume of buffer and centrifuge at $300 \times g$ for 10 minutes at $4-8^\circ\text{C}$. Pipette off supernatant completely.
7. Resuspend cell pellet in buffer:
Depletion with LD Column: 500 μL for up to 1.25×10^8 cells
Depletion with autoMACS Separator: 500 μL for up to 1×10^8 cells
▲ **Note:** For larger cell numbers, scale up buffer volume accordingly.
8. Proceed to magnetic separation (2.3).



2.3 Magnetic separation: Depletion of non-T cells

Depletion with LD Column

1. Place LD Column in the magnetic field of a suitable MACS Separator (see "Column data sheets").
2. Prepare column by rinsing with 2 mL of buffer.
3. Apply cell suspension onto the column.
4. Collect unlabeled cells which pass through and wash column with 2×1 mL of buffer. Perform washing steps by adding buffer successively once the column reservoir is empty. Collect total effluent. This contains the unlabeled pre-enriched T cell fraction.
5. Proceed to 2.4 for the isolation of TCR γ/δ^+ T cells.

Depletion with the autoMACS™ Separator

▲ Refer to the "autoMACS™ User Manual" for instructions on how to use the autoMACS Separator.

1. Prepare and prime the autoMACS Separator.
2. Place the tube containing the magnetically labeled cells in the autoMACS Separator. Choose separation program "Depl05".
3. Collect the unlabeled fraction (outlet port "neg1"). This is the pre-enriched T cell fraction.
4. Proceed to 2.4 for the isolation of TCR γ/δ^+ T cells.



2.4 Magnetic labeling of TCR γ/δ^+ T cells

▲ Volumes for magnetic labeling given below are for an initial starting cell number of up to 10^8 leukocytes. For larger initial cell numbers, scale up volumes accordingly.

1. Centrifuge the cells at $300 \times g$ for 10 minutes. Pipette off supernatant completely.
2. Resuspend cell pellet in 450 μL of buffer.

3. Add 50 μL of **Anti-Biotin MicroBeads**.
4. Mix well and incubate for 15 minutes at 4–8 °C.
5. Wash cells by adding 10–20 $\times$ labeling volume of buffer and centrifuge at 300 $\times$ g for 10 minutes. Pipette off supernatant completely.
6. Resuspend up to 10^8 cells in 500 μL of buffer.
7. Proceed to magnetic separation (2.5).



2.5 Magnetic separation: Positive selection of TCR γ/δ^+ T cells

Positive selection with MS Columns

1. Place MS Column in the magnetic field of a suitable MACS Separator (see "Column data sheets").
 2. Prepare column by rinsing with 500 μL of buffer.
 3. Apply cell suspension onto the column.
 4. Wash column with 3 $\times$ 500 μL of buffer. Perform washing steps by adding buffer three times once the column reservoir is empty.
 5. Remove column from the separator and place it on a suitable collection tube.
 6. Pipette 1 mL of buffer onto the column. Immediately flush out the fraction with magnetically labeled TCR γ/δ^+ T cells by firmly applying the plunger supplied with the column.
- iii) Repeat the magnetic separation procedure as described in steps 1 to 6 by using a new column.


Positive selection with the autoMACS™ Separator

▲ Refer to the "autoMACS™ User Manual" for instructions on how to use the autoMACS Separator.

1. Prepare and prime the autoMACS Separator.
2. Place the tube containing the magnetically labeled cells in the autoMACS Separator. Choose separation program "Posseld2".
3. Collect the positive fraction (outlet port "pos2"). This is the enriched TCR γ/δ^+ T cell fraction.

3. Example of a separation using the TCR γ/δ^+ T Cell Isolation Kit

TCR γ/δ^+ T cells were isolated from single-cell suspensions of mouse spleen (A-C), lymph nodes (D), and thymus (E) using the TCR γ/δ^+ T Cell Isolation Kit, an LD and two MS Columns, a MidiMACS™ and a MiniMACS™ Separator. The cells were fluorescently stained with CD3-FITC, Anti-TCR γ/δ -Biotin and Anti-Biotin-APC (# 130-090-856). Cell debris and dead cells were excluded from the analysis based on scatter signals and PI fluorescence.

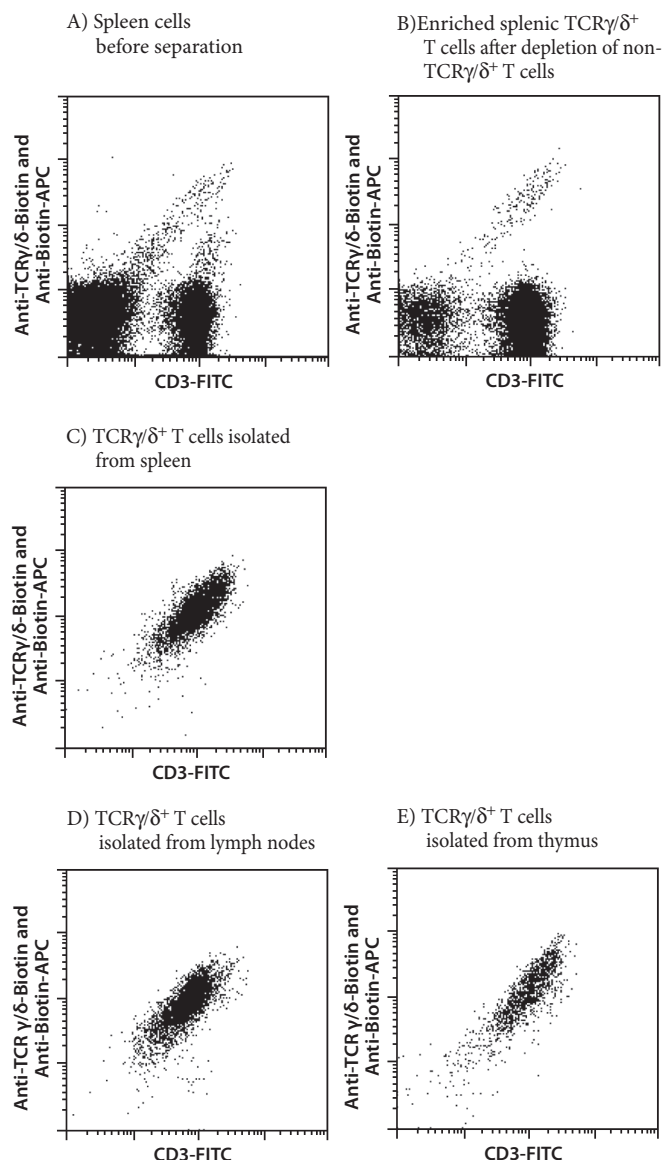

4. Reference

1. Carding, SR; Egan, PJ (2002) Gammadelta T cells: functional plasticity and heterogeneity. Nat. Rev. Immunol. 5: 336–345.

Refer to www.miltenyibiotec.com for all data sheets and protocols. Miltenyi Biotec provides technical support worldwide. Visit www.miltenyibiotec.com/local to find your nearest Miltenyi Biotec contact.

Warnings

Reagents contain sodium azide. Under acidic conditions sodium azide yields hydrazoic acid, which is extremely toxic. Azide compounds should be diluted with running water before discarding. These precautions are recommended to avoid deposits in plumbing where explosive conditions may develop.

Legal notices

Limited product warranty

Miltenyi Biotec B.V. & Co. KG and/or its affiliate(s) warrant this product to be free from material defects in workmanship and materials and to conform substantially with Miltenyi Biotec's published specifications for the product at the time of order, under normal use and conditions in accordance with its applicable documentation, for a period beginning on the date of delivery of the product by Miltenyi Biotec or its authorized distributor and ending on the expiration date of the product's applicable shelf life stated on the product label, packaging or documentation (as applicable) or, in the absence thereof, ONE (1) YEAR from date of delivery ("Product Warranty"). Miltenyi Biotec's Product Warranty is provided subject to the warranty terms as set forth in Miltenyi Biotec's General Terms and Conditions for the Sale of Products and Services available on Miltenyi Biotec's website at www.miltenyibiotec.com, as in effect at the time of order ("Product Warranty"). Additional terms may apply. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING IF A PRODUCT IS SUITABLE FOR CUSTOMER'S PARTICULAR PURPOSE AND APPLICATION METHODS.

Technical information

The technical information, data, protocols, and other statements provided by Miltenyi Biotec in this document are based on information, tests, or experience which Miltenyi Biotec believes to be reliable, but the accuracy or completeness of such information is not guaranteed. Such technical information and data are intended for persons with knowledge and technical skills sufficient to assess and apply their own informed judgment to the information. Miltenyi Biotec shall not be liable for any technical or editorial errors or omissions contained herein.

All information and specifications are subject to change without prior notice. Please contact Miltenyi Biotec Technical Support or visit www.miltenyibiotec.com for the most up-to-date information on Miltenyi Biotec products.

Licenses

This product and/or its use may be covered by one or more pending or issued patents and/or may have certain limitations. Certain uses may be excluded by separate terms and conditions. Please contact your local Miltenyi Biotec representative or visit Miltenyi Biotec's website at www.miltenyibiotec.com for more information.

The purchase of this product conveys to the customer the non-transferable right to use the purchased amount of the product in research conducted by the customer (whether the customer is an academic or for-profit entity). This product may not be further sold. Additional terms and conditions (including the terms of a Limited Use Label License) may apply.

CUSTOMER'S USE OF THIS PRODUCT MAY REQUIRE ADDITIONAL LICENSES DEPENDING ON THE SPECIFIC APPLICATION. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING FOR ITSELF WHETHER IT HAS ALL APPROPRIATE LICENSES IN PLACE. Miltenyi Biotec provides no warranty that customer's use of this product does not and will not infringe intellectual property rights owned by a third party. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

Trademarks

autoMACS, MACS, MidiMACS, the Miltenyi Biotec logo, MiniMACS, OctoMACS, QuadroMACS, SuperMACS and VarioMACS are registered trademarks or trademarks of Miltenyi Biotec and/or its affiliates in various countries worldwide.

Copyright © 2020 Miltenyi Biotec and/or its affiliates. All rights reserved.