

Index

1. Description
 - 1.1 Principle of MACS® separation
 - 1.2 Background and product applications
 - 1.3 Reagent and instrument requirements
2. Protocol
 - 2.1 Sample preparation
 - 2.2 Magnetic labeling
 - 2.3 Magnetic separation
3. Example of a separation using Anti-DC (OX62) MicroBeads
4. Appendix
5. References

1. Description

Components	2 mL Anti-DC (OX62) MicroBeads, rat: MicroBeads conjugated to monoclonal mouse anti-DC (OX62) antibodies (isotype: mouse IgG1; clone: OX62).
Size	For 10 ⁹ total cells, up to 100 separations.
Product format	The product is supplied as suspension containing 0.1% gelatine and 0.05% sodium azide.
Storage	Store protected from light at 4–8 °C. Do not freeze. The expiration date is indicated on the vial label.

1.1 Principle of MACS® separation

First the OX62⁺ cells are magnetically labeled with Anti-DC (OX62) MicroBeads. Then the cell suspension is loaded onto a MACS® Column which is placed in the magnetic field of a MACS Separator. The magnetically labeled OX62⁺ cells are retained on the column. The unlabeled cells run through and this cell fraction is depleted of OX62⁺ cells. After removal of the column from the magnetic field, the magnetically retained OX62⁺ cells can be eluted as the positively selected cell fraction.

1.2 Background and product applications

Anti-DC (OX62) MicroBeads are developed for the positive selection or depletion of OX62⁺ dendritic cells (DC) from blood, lymphoid and non-lymphoid tissue. The monoclonal antibody clone OX62 recognizes the integrin α_{E2} subunit¹ (molecular mass 150 kDa)² expressed on dendritic cells in the rat. Notably, OX62 antigen is found on both CD4⁺ OX41⁺ and CD4⁻ OX41⁻ dendritic cells.^{3,4,5} The OX62 antigen was also shown to be expressed on MHC II⁺ TCR γ / δ ⁺ T cells¹, and on a subset of folliculo-stellate like (S100⁺) cells in the anterior pituitary^{6,7}. OX62 is not detectable on Langerhans cells.¹

Examples of applications

- Positive selection or depletion of dendritic cells from single-cell suspensions of lymphoid or non-lymphoid tissue, or peripheral blood.⁸
- ### 1.3 Reagent and instrument requirements
- Buffer (degassed): Prepare a solution containing PBS (phosphate buffered saline) pH 7.2, 0.5% BSA and 2 mM EDTA by diluting MACS BSA Stock Solution (#130-091-376) 1:20 in autoMACS™ Rinsing Solution (# 130-091-222). Keep buffer cold (4–8 °C).

▲ **Note:** EDTA can be replaced by other supplements such as anticoagulant citrate dextrose formula-A (ACD-A) or citrate phosphate dextrose (CPD). BSA can be replaced by other proteins such as rat serum albumin, rat serum or fetal calf serum. Buffers or media containing Ca²⁺ or Mg²⁺ are not recommended for use.
 - MACS Columns and MACS Separators: OX62⁺ cells can be enriched by using MS, LS or XS Columns (positive selection). Anti-DC (OX62) MicroBeads can be used for depletion of OX62⁺ cells on LD, CS or D Columns. Cells which strongly express the OX62 antigen can also be depleted using MS, LS or XS Columns. Positive selection or depletion can also be performed by using the autoMACS Separator.

Column	max. number of labeled cells	max. number of total cells	Separator
Positive selection			
MS	10 ⁷	2×10 ⁸	MiniMACS, OctoMACS, VarioMACS, SuperMACS
LS	10 ⁸	2×10 ⁹	MidiMACS, QuadroMACS, VarioMACS, SuperMACS
XS	10 ⁹	2×10 ¹⁰	SuperMACS
Depletion			
LD	10 ⁸	5×10 ⁸	MidiMACS, QuadroMACS, VarioMACS, SuperMACS
CS	2×10 ⁸		VarioMACS, SuperMACS
D	10 ⁹		SuperMACS
Positive selection or depletion			
autoMACS	2×10 ⁸	4×10 ⁹	autoMACS

▲ **Note:** Column adapters are required to insert certain columns into VarioMACS™ Separator or SuperMACS™ Separator. For details, see MACS Separator data sheets.

- Collagenase D: 2 mg/mL Collagenase D (>0.15 U/ mg, e.g. from Roche, Germany) in RPMI/ 1% FCS.
- (Optional) Nycodenz®-PBS solution. Nycodenz® (from Nycomed As, Oslo, Norway) solubilized in PBS buffer at a final concentration of 14.5% (w/v).
- (Optional) Fluorochrome-conjugated Anti-DC (OX62) antibody for flow-cytometric analysis.

- (Optional) PI (propidium iodide) or 7-AAD for flow-cytometric exclusion of dead cells.
- (Optional) Pre-Separation Filters (# 130-041-407) to remove cell clumps.

2. Protocol

2.1 Sample preparation

Prepare a single-cell suspension from lymphoid organs, non-lymphoid tissue or peripheral blood using standard methods (see "General Protocols" in the User Manuals or visit www.miltenyibiotec.com/protocols).

▲ **Note:** Dead cells may bind non-specifically to MACS MicroBeads. In case of high numbers of dead cells, removal of dead cells by density gradient centrifugation or the Dead Cell Removal Kit (# 130-090-101) is recommended.

Preparation of rat spleens with Collagenase D

To increase the recovery and final purity of dendritic cells, it is recommended to prepare spleen cell suspensions by Collagenase D treatment:

1. Place up to 2 isolated spleens in a 6 cm petri-dish with sufficient Collagenase D to completely cover the bottom of the dish (about 5 mL).
2. Inject spleens with 500 μ L of Collagenase D per spleen using a 1 mL syringe and a 25G needle and cut the tissue in smaller pieces using a pair of scissors and a pair of forceps.
3. Incubate spleen pieces for 25 minutes at 37 °C in Collagenase D. Add 10 mM EDTA and incubate a further 5 minutes.
4. Pass the digested material through a steel mesh using a plunger of a syringe and wash the petri-dish and the mesh with buffer.
5. Pass all cells through a Pre-Separation Filters (# 130-041-407) to remove clumps and collect cells in a 50 mL tube. Add buffer to a final volume of about 50 mL.
6. Wash cells by centrifugation at 200 \times g for 10 minutes. Count cells.
7. Proceed to Nycodenz gradient centrifugation or magnetic labeling (2.2)

(Optional) Pre-enrichment of rat dendritic cells by Nycodenz® gradient centrifugation

For highly pure dendritic cells, OX62⁺ CD3⁺ TCR γ/δ ⁺ T cells might be depleted prior to OX62 separation by Nycodenz® gradient centrifugation.

1. After centrifugation remove supernatant completely and resuspend cell pellet in 1 mL of buffer per 1–2 \times 10⁸ total cells. For fewer cells, use same volume.
2. Layer cell suspension on equal volume of 14.5% (w/v) Nycodenz-PBS solution (1.079 g/mL) and centrifuge for 13 minutes at 1,800 \times g and 4 °C.
3. Recover the low density cells at the top of the low density solution. Add buffer to a final volume of 50 mL and wash cells by centrifugation for 10 minutes at 200 \times g and 4 °C. Count cells.



2.2 Magnetic labeling

▲ Work fast, keep cells cold, and use pre-cooled solutions. This will prevent capping of antibodies on the cell surface and non-specific cell labeling.

▲ Volumes for magnetic labeling given below are for up to 10⁷ total cells. When working with fewer than 10⁷ cells, use the same volumes as indicated. When working with higher cell numbers, scale up all reagent volumes and total volumes accordingly (e.g. for 2 \times 10⁷ total cells, use twice the volume of all indicated reagent volumes and total volumes).

▲ For optimal performance it is important to obtain a single-cell suspension before magnetic separation. Pass cells through 30 μ m nylon mesh (Pre-Separation Filters # 130-041-407) to remove cell clumps which may clog the column.

1. Determine cell number.
2. Centrifuge cell suspension at 300 \times g for 10 minutes. Pipette off supernatant completely.
3. Resuspend cell pellet in 80 μ L of buffer per 10⁷ total cells.
4. Add 20 μ L of Anti-Rat DC (OX62) MicroBeads per 10⁷ total cells.
5. Mix well and incubate for 15 minutes at 4–8 °C.
 - ▲ **Note:** Working on ice may require increased incubation times. Higher temperatures and/or longer incubation times lead to non-specific cell labeling.
6. (Optional) Add staining antibodies and incubate for 5 minutes at 4–8 °C.
7. Wash cells by adding 1–2 mL of buffer per 10⁷ cells and centrifuge at 300 \times g for 10 minutes. Pipette off supernatant completely.
8. Resuspend up to 10⁸ cells in 500 μ L of buffer.
 - ▲ **Note:** For higher cell numbers, scale up buffer volume accordingly.
 - ▲ **Note:** For depletion with LD Columns, resuspend up to 1.25 \times 10⁸ cells in 500 μ L of buffer.
9. Proceed to magnetic separation (2.3).



2.3 Magnetic separation

▲ Choose an appropriate MACS Column and MACS Separator according to the number of total cells and the number of OX62⁺ cells (see table in section 1.3).

Magnetic separation with MS or LS Columns

1. Place column in the magnetic field of a suitable MACS Separator (see "Column data sheets").
2. Prepare column by rinsing with appropriate amount of buffer:

MS: 500 μ L	LS: 3 mL.
-----------------	-----------
3. Apply cell suspension onto the column.
4. Collect unlabeled cells which pass through and wash column with appropriate amount of buffer. Perform washing steps by adding buffer three times, each time once the column reservoir is empty.

MS: 3 \times 500 μ L	LS: 3 \times 3 mL.
----------------------------	----------------------

 Collect total effluent. This is the unlabeled cell fraction.

- Remove column from the separator and place it on a suitable collection tube.
- Pipette appropriate amount of buffer onto the column. Immediately flush out fraction with the magnetically labeled cells by firmly applying the plunger supplied with the column.

MS: 1 mL LS: 5 mL.

▲ **Note:** To increase the purity of the magnetically labeled fraction, it can be passed over a new, freshly prepared column.

Magnetic separation with XS Columns

For instructions on the column assembly and the separation, refer to the "XS Column data sheet".

Depletion with LD Columns

- Place LD Column in the magnetic field of a suitable MACS Separator (see "LD Column data sheet").
- Prepare column by rinsing with 2 mL of buffer.
- Apply cell suspension onto the column.
- Collect unlabeled cells which pass through and wash column with 2×1 mL of buffer. Collect total effluent. This is the unlabeled cell fraction.

Depletion with CS Columns

- Assemble CS Column and place it in the magnetic field of a suitable MACS Separator (see "CS Column data sheet").
- Prepare column by filling and rinsing with 60 mL of buffer. Attach a 22G flow resistor to the 3-way-stopcock of the assembled column (see "CS Column data sheet").
- Apply cell suspension onto the column.
- Collect unlabeled cells which pass through and wash column with 30 mL buffer from the top. Collect total effluent. This is the unlabeled cell fraction.

Depletion with D Columns

For instructions on column assembly and separation, refer to the "D Column data sheet".

Magnetic separation with the autoMACS™ Separator

▲ Refer to the "autoMACS™ User Manual" for instructions on how to use the autoMACS Separator.

- Prepare and prime autoMACS Separator.
- Place tube containing the magnetically labeled cells in the autoMACS Separator. For a standard separation, choose following separation programs:

Positive selection: "Possel"

Depletion: "Depletes"


▲ **Note:** Program choice depends on the isolation strategy, the strength of magnetic labeling and the frequency of magnetically labeled cells. For details see autoMACS User Manual: "autoMACS Cell Separation Programs".

- When using the program "Possel", collect positive fraction (outlet port "pos1"). This is the purified OX62⁺ cell fraction.


When using the program "Depletes", collect unlabeled fraction (outlet port "neg1"). This is the OX62⁻ cell fraction.

3. Example of a separation using Anti-DC (OX62) MicroBeads


Dendritic cells were pre-enriched by Nycodenz gradient centrifugation from rat spleen cell suspension and were then isolated using Anti-DC (OX62) MicroBeads and a MiniMACS™ Separator with an MS Column. For flow cytometric analysis, cells were stained for OX62 and MHC Class II (OX6). Cell debris and dead cells were excluded from the analysis based on scatter signals and PI fluorescence.

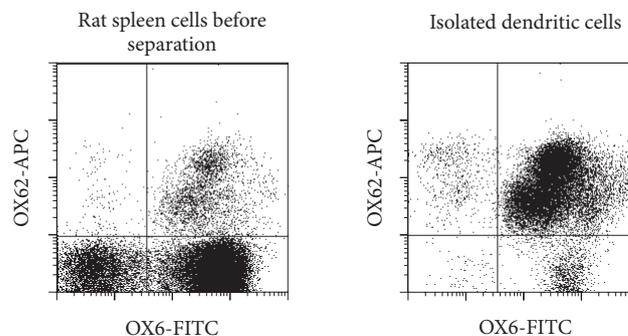

To evaluate the depletion of OX62⁺CD3⁺TCR γ/δ T cells from rat spleen by Nycodenz-preparation followed by MACS Technology, a sample of the isolated dendritic cells was stained with CD3-PE.

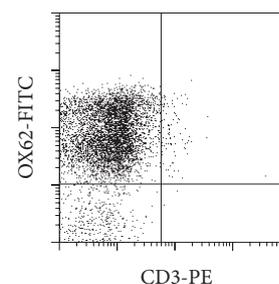

To distinguish between the two major OX62⁺ dendritic cells subsets a sample of the isolated dendritic cells was additionally stained for CD4 (OX35). Note that OX62^{bright} dendritic cells are CD4⁻, whereas OX62^{dim} cells are CD4⁺.

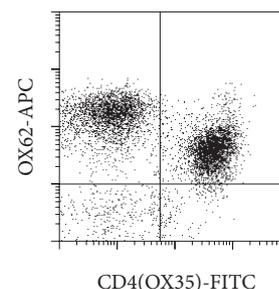

5. References

1. Brenan, M; Rees, JG (1997) Sequence analysis of rat integrin α_{E1} and α_{E2} subunits. tissue expression reveals phenotypic similarities between intraepithelial lymphocytes and dendritic cells in lymph. *Eur. J. Immunol.* 27: 3070-3079. [957]
2. Brenan, M; Puklavec, M (1992) The MRC OX-62 Antigen: A Useful Marker in the Purification of Rat Veiled Cells with the Biochemical Properties of an Integrin. *J. Exp. Med.* 175: 1457-1465. [14]
3. Huang, F-P; Platt, N; Wykes, M; Major, JR; Powell, TJ; Jenkins, CD; MacPherson, GG (2000) A Discrete Subpopulation of Dendritic Cells Transports Apoptotic Intestinal Epithelial Cells to T Cell Areas of Mesenteric Lymph Nodes. *J. Exp. Med.* 191: 435-443. [953]
4. Trinité, B; Voisine, C; Yagita, H; Josien, R (2000) A Subset of Cytolytic Dendritic Cells in Rat. *J. Immunol.* 165: 4202-4208. [955]
5. Liu, L; Zhang, M; Jenkins, C; MacPherson, GG (1998) Dendritic Cell Heterogeneity In Vivo: Two Functionally Different Dendritic Cell Populations in Rat Intestinal Lymph Can Be Distinguished by CD4 Expression. *J. Immunol.* 161: 1146-1155.
6. Allaerts, W; Jeucken, PHM; Debets, R; Hoefakker, S; Claassen, E; Drexhage, HA (1997) Heterogeneity of Pituitary Folliculo-Stellate Cells: Implications for Interleukin-6 Production and Accessory Function *in vitro*. *J. Neuroendocrinol.* 9: 43-53. [956]
7. Allaerts, W; Salomon, B; Leenen, PJM; van Wijngaardt, S; Jeucken, PHM; Ruuls, S; Klatzmann, D; Drexhage, HA (1997) A population of interstitial cells in the anterior pituitary with a hematopoietic origin and a rapid turnover: a relationship with folliculo-stellate cells? *J. Neuroimmunol.* 78: 184-197.
8. Kobayashi H, Miura S, Nagata H, Tsuzuki Y, Hokari R, Ogino T, Watanabe C, Azuma T, Ishii H. (2004) *In situ* demonstration of dendritic cell migration from rat intestine to mesenteric lymph nodes: relationships to maturation and role of chemokines. *J. Leukoc. Biol.* 75(3): 434-42. [3995]

Refer to www.miltenyibiotec.com for all data sheets and protocols. Miltenyi Biotec provides technical support worldwide. Visit www.miltenyibiotec.com/local to find your nearest Miltenyi Biotec contact.

Warnings

Reagents contain sodium azide. Under acidic conditions sodium azide yields hydrazoic acid, which is extremely toxic. Azide compounds should be diluted with running water before discarding. These precautions are recommended to avoid deposits in plumbing where explosive conditions may develop.

Legal notices

Limited product warranty

Miltenyi Biotec B.V. & Co. KG and/or its affiliate(s) warrant this product to be free from material defects in workmanship and materials and to conform substantially with Miltenyi Biotec's published specifications for the product at the time of order, under normal use and conditions in accordance with its applicable documentation, for a period beginning on the date of delivery of the product by Miltenyi Biotec or its authorized distributor and ending on the expiration date of the product's applicable shelf life stated on the product label, packaging or documentation (as applicable) or, in the absence thereof, ONE (1) YEAR from date of delivery ("Product Warranty"). Miltenyi Biotec's Product Warranty is provided subject to the warranty terms as set forth in Miltenyi Biotec's General Terms and Conditions for the Sale of Products and Services available on Miltenyi Biotec's website at www.miltenyibiotec.com, as in effect at the time of order ("Product Warranty"). Additional terms may apply. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING IF A PRODUCT IS SUITABLE FOR CUSTOMER'S PARTICULAR PURPOSE AND APPLICATION METHODS.

Technical information

The technical information, data, protocols, and other statements provided by Miltenyi Biotec in this document are based on information, tests, or experience which Miltenyi Biotec believes to be reliable, but the accuracy or completeness of such information is not guaranteed. Such technical information and data are intended for persons with knowledge and technical skills sufficient to assess and apply their own informed judgment to the information. Miltenyi Biotec shall not be liable for any technical or editorial errors or omissions contained herein.

All information and specifications are subject to change without prior notice. Please contact Miltenyi Biotec Technical Support or visit www.miltenyibiotec.com for the most up-to-date information on Miltenyi Biotec products.

Licenses

This product and/or its use may be covered by one or more pending or issued patents and/or may have certain limitations. Certain uses may be excluded by separate terms and conditions. Please contact your local Miltenyi Biotec representative or visit Miltenyi Biotec's website at www.miltenyibiotec.com for more information.

The purchase of this product conveys to the customer the non-transferable right to use the purchased amount of the product in research conducted by the customer (whether the customer is an academic or for-profit entity). This product may not be further sold. Additional terms and conditions (including the terms of a Limited Use Label License) may apply.

CUSTOMER'S USE OF THIS PRODUCT MAY REQUIRE ADDITIONAL LICENSES DEPENDING ON THE SPECIFIC APPLICATION. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING FOR ITSELF WHETHER IT HAS ALL APPROPRIATE LICENSES IN PLACE. Miltenyi Biotec provides no warranty that customer's use of this product does not and will not infringe intellectual property rights owned by a third party. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

Trademarks

autoMACS, MACS, MidiMACS, the Miltenyi Biotec logo, MiniMACS, OctoMACS, QuadroMACS, SuperMACS, and VarioMACS are registered trademarks or trademarks of Miltenyi Biotec and/or its affiliates in various countries worldwide.

Copyright © 2020 Miltenyi Biotec and/or its affiliates. All rights reserved.