

Contents

1. Description
 - 1.1 Principle of the NK Cell Activation/Expansion Kit, human
 - 1.2 Background information
 - 1.3 Applications
 - 1.4 Reagent and instrument requirements
2. Protocol
 - 2.1 Loading of MACS Anti-Biotin MACSiBead Particles
 - 2.2 Magnetic separation of NK cells using the NK Cell Isolation Kit, human
 - 2.3 NK cell activation and expansion protocol
 - 2.4 Immunofluorescent staining
3. Examples of NK cell activation and expansion using the NK Cell Activation/Expansion Kit, human

1. Description

This product is for research use only.

Components	<p>2 mL MACS® Anti-Biotin MACSiBead™ Particles, cell culture grade, corresponding to 4×10⁸ MACSiBead Particles; MACSiBead Particles conjugated to monoclonal anti-biotin antibodies.</p> <p>0.4 mL CD335 (NKp46)-Biotin, human – functional grade (100 µg/mL).</p> <p>0.4 mL CD2-Biotin, human – functional grade (100 µg/mL).</p>
Product format	All reagents are supplied in azide-free buffer, MACS Anti-Biotin MACSiBead Particles contain stabilizer.
Storage	Store protected from light at 2–8 °C. Do not freeze. The expiration date is indicated on the vial label.

1.1 Principle of the NK Cell Activation/Expansion Kit, human

The NK Cell Activation/Expansion Kit is designed to activate and expand human NK cells. The kit consists of MACS Anti-Biotin MACSiBead Particles and biotinylated antibodies against human CD335 (NKp46) and CD2. MACS Anti-Biotin MACSiBead Particles loaded with biotinylated antibodies are used to activate and expand resting NK cells purified from human blood or peripheral blood mononuclear cells (PBMCs).

1.2 Background information

In a first step the MACS Anti-Biotin MACSiBead Particles are loaded with biotinylated antibodies. Best activation is achieved by using equal amounts of the provided biotinylated antibodies against CD335 (NKp46) and CD2.

▲ **Note:** Other combinations of biotinylated antibodies may be experimentally tested for their suitability, if required.

Loaded MACS Anti-Biotin MACSiBead Particles are subsequently used for the expansion of NK cells. Best activation and expansion of NK cells is accomplished by using one loaded MACS Anti-Biotin MACSiBead Particle per two cells (bead-to-cell ratio 1:2). The cells are cultured for further expansion.

NK cells, activated by using MACS Anti-Biotin MACSiBead Particles, can be used for any downstream processing such as cytokine analysis, cytolytic activity, gene expression, or functional studies.

MACS Anti-Biotin MACSiBead Particles show no autofluorescence and normally do not need to be removed prior to flow cytometric analysis.

1.3 Applications

- Activation and expansion of resting NK cells from PBMCs or after isolation with the NK Cell Isolation Kit, human (# 130-092-657) or CD56 MicroBeads, human (# 130-050-401).

1.4 Reagent and instrument requirements

- Buffer: Phosphate-buffered saline (PBS) pH 7.2, supplemented with 0.5% human serum albumin (HSA) and 2 mM EDTA. Keep buffer cold (2–8 °C). Degas buffer before use, as air bubbles could block the column.

▲ **Note:** HSA can be replaced by other proteins such as bovine serum albumin, fetal calf serum or human AB serum. Buffers or media containing Ca²⁺ or Mg²⁺ are not recommended for use.
- NK Cell Isolation Kit, human (# 130-092-657) or CD56 MicroBeads, human (# 130-050-401).
- MACS Columns and MACS Separators: Choose the appropriate MACS Separator and MACS Columns according to the number of labeled cells and to the number of total cells.

Column	Max. number of labeled cells	Max. number of total cells	Separator
Positive selection			
MS	10 ⁷	2×10 ⁸	MiniMACS, OctoMACS, SuperMACS II
LS	10 ⁸	2×10 ⁹	MidiMACS, QuadroMACS, SuperMACS II

▲ **Note:** Column adapters are required to insert certain columns into SuperMACS™ II Separators. For details refer to the respective MACS Separator data sheet.

- Medium: NK MACS Medium, human (# 130-114-429) supplemented with 5% AB serum and 500 IU/mL human Interleukin 2 (IL-2).

- Human IL-2, e.g., Human IL-2 IS, premium grade (# 130-097-748).
- (Optional) MACS GMP Cell Expansion Bags (# 170-076-403) or flat bottom cell culture plates with lids.
- Humidified incubator.
- MACSmix™ Tube Rotator (# 130-090-753) for loading of MACSiBead Particles.
- (Optional) Fluorochrome-conjugated antibodies for flow cytometric analysis. For more information about antibodies refer to www.miltenyibiotec.com/antibodies.
- (Optional) Propidium Iodide Solution (# 130-093-233) or 7-AAD Staining Solution (# 130-111-568) for flow cytometric exclusion of dead cells.

2. Protocol

▲ All steps in the protocol have to be performed under sterile conditions.

2.1 Loading of MACS Anti-Biotin MACSiBead Particles

▲ Resuspend MACS Anti-Biotin MACSiBead Particles thoroughly by vortexing before use, to obtain a homogenous suspension.

▲ MACS Anti-Biotin MACSiBead Particles are supplied without preservative. Remove aliquots under aseptic conditions.

▲ It is recommended to load MACS Anti-Biotin MACSiBead Particles in batches of 1×10^8 MACS Anti-Biotin MACSiBead Particles. Loaded MACS Anti-Biotin MACSiBead Particles are stable for up to two months when stored at 2–8 °C.

1. Pipette 100 µL of CD335 (NKp46)-Biotin and 100 µL CD2-Biotin into sealable 2 mL tube and mix well.

▲ **Note:** This antibody combination is optimized for achieving maximum NK cell activation and expansion.

2. Resuspend MACS Anti-Biotin MACSiBead Particles thoroughly by vortexing.
3. Remove 500 µL MACS Anti-Biotin MACSiBead Particles (1×10^8 MACS Anti-Biotin MACSiBead Particles) and add to antibody mix.
4. Add 300 µL buffer to adjust to a total volume of 1 mL.

▲ **Note:** MACS Anti-Biotin MACSiBead Particles can be loaded in a flexible manner with biotinylated antibodies or ligands other than those supplied. If desired, add other biotinylated antibodies or ligands at appropriate concentrations and adjust with buffer to a total volume of 1 mL, accordingly.
5. Incubate for 2 hours at 2–8 °C under constant, gentle rotation by using the MACSmix Tube Rotator at approximately 12 rpm (slowest permanent run program).
6. The loaded MACS Anti-Biotin MACSiBead Particles (1×10^8 MACS Anti-Biotin MACSiBead Particles/mL) are now ready to use. Do not remove the loaded MACS Anti-Biotin MACSiBead Particles from the antibody mix. Store at 2–8 °C for up to two months.



2.2 Magnetic separation of NK cells using the NK Cell Isolation Kit, human

▲ Isolate NK cells according to the data sheet of the NK Cell Isolation Kit, human. The MACS Columns and MACS Separators that can be used in this step are listed above (see 1.4).

2.3 NK cell activation and expansion protocol

This NK cell activation and expansion protocol is optimized for NK cells that have been purified using the NK Cell Isolation Kit, human using one loaded MACS Anti-Biotin MACSiBead Particle per two NK cells (bead-to-cell ratio 1:2).

▲ **Note:** Ratios other than 1:2 of loaded MACS Anti-Biotin MACSiBead Particles per cell may be required for other applications.

▲ Volumes for activation given below are for 10^6 purified NK cells or 10^6 PBMCs. When working with higher cell numbers, scale up all reagent volumes and total volumes accordingly (e.g. for 2×10^6 total cells, use twice the volume of all indicated reagent volumes and total volumes).

1. Resuspend loaded MACS Anti-Biotin MACSiBead Particles thoroughly and transfer 5 µL (5×10^5 loaded MACS Anti-Biotin MACSiBead Particles) per 10^6 NK cells to a suitable tube.

▲ **Note:** If unloaded MACS Anti-Biotin MACSiBead Particles shall be used for negative control experiments, replace loaded MACS Anti-Biotin MACSiBead Particles by adding 5×10^5 of unloaded MACS Anti-Biotin MACSiBead Particles per 10^6 NK cells.

2. Add 100 µL culture medium to the loaded MACS Anti-Biotin MACSiBead Particles and centrifuge at 300×g for 5 minutes.
3. Aspirate supernatant and resuspend loaded MACS Anti-Biotin MACSiBead Particles in 50 µL of fresh culture medium.
4. Resuspend PBMCs or purified NK cells at a density of 10^6 cells per 950 µL of culture medium (NK MACS Medium supplemented with 5% AB serum and 500 IU/mL IL-2).
5. Add the prepared MACS Anti-Biotin MACSiBead Particles from step 3 to the 950 µL of cell suspension and mix well.
6. Add the mixture to a suitable cell culture vessel at a density of 10^6 cells per mL, for example, in the wells of a 24-well cell culture plate.
7. Incubate at 37 °C and 5% CO₂.

▲ **Note:** Inspect cultures daily, and add fresh medium if required.

▲ **Note:** For NK cell expansion the addition of culture medium is required regularly. In the following a guideline for the stimulation and expansion of NK cells is described.

8. At day 6, gently pipette cell suspension up and down to break up clumps.
9. Determine cell number and dilute to $1\text{--}1.5 \times 10^6$ cells per mL by adding fresh culture medium (NK MACS Medium supplemented with 5% AB serum and 500 IU/mL IL-2). Transfer to a fresh culture vessel of appropriate size.

▲ **Note:** NK cell expansion is donor-dependent. Best results are achieved when NK cells are maintained in a cell density of $1\text{--}1.5 \times 10^6$ NK cells per mL. Depending on the expansion rate it might be necessary to split culture daily.

2.4 Immunofluorescent staining

▲ Volumes for fluorescent labeling given below are for 10^6 total cells. When working with fewer than 10^6 cells and up to 10^7 cells, use the same volumes as indicated.

▲ MACS Anti-Biotin MACSiBead Particles show no autofluorescence and do not need to be removed prior to flow cytometric analysis.

▲ Scatter properties of cells may be altered due to strong interaction between cells and MACS Anti-Biotin MACSiBead Particles.

1. Resuspend cells to break up cell clumps.
2. Wash cells by adding 1–2 mL of buffer per 10^6 cells and centrifuge at $300\times g$ for 10 minutes. Aspirate supernatant completely.
3. Add 10 μ L of each staining antibody, e.g., CD3 Antibody, anti-human, FITC and CD56 Antibody, anti-human, PE to 10^6 cells resuspended with buffer to a total volume of 110 μ L.
4. Mix well and incubate for 10 minutes in the dark in the refrigerator (2–8 °C).


▲ **Note:** Working on ice requires increased incubation time. Higher temperatures and/or longer incubation times lead to non-specific cell labeling.

5. Wash cells by adding 1–2 mL of buffer per 10^6 cells and centrifuge at $300\times g$ for 10 minutes. Aspirate supernatant completely.
6. Resuspend cell pellet in a suitable amount of buffer for analysis by flow cytometry or fluorescence microscopy.

3. Example of NK cell activation and expansion using the NK Cell Activation/Expansion Kit, human

A) Cytotoxicity of expanded NK cells


MACS Anti-Biotin MACSiBead Particles were loaded with CD335 (NKp46) and CD2 antibodies. NK cells were activated and expanded from PBMCs using one loaded MACS Anti-Biotin MACSiBead Particle per viable cell. Cells were cultured in NK MACS Medium supplemented with 5% AB serum and 500 IU/mL IL-2. After two weeks, NK cells were incubated with K562 target cells in different ratios. In addition, NK cells incubated in NK MACS Medium supplemented with 5% AB serum and 500 IU/mL IL-2 alone were analyzed. Specific target cell lysis by NK cells was determined after 4 hours.

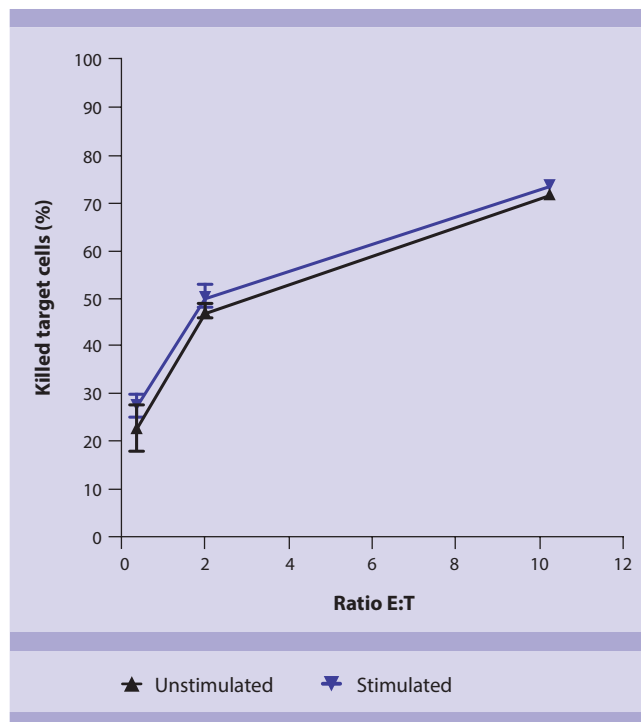

B) NK cell expansion rates

MACS Anti-Biotin MACSiBead Particles were loaded with CD335 (NKp46) and CD2 antibodies. NK cells were isolated using the NK Cell Isolation Kit, human and expanded using one loaded MACS Anti-Biotin MACSiBead Particle per two NK cells. Cells were cultured in NK MACS Medium supplemented with 5% AB serum and 500 IU/mL IL-2 at an initial density of 10^6 NK cells per mL. Cells were expanded for 18 days.

For comparison NK cells were cultured in medium supplemented with 5% AB serum and 500 IU/mL IL-2 alone.

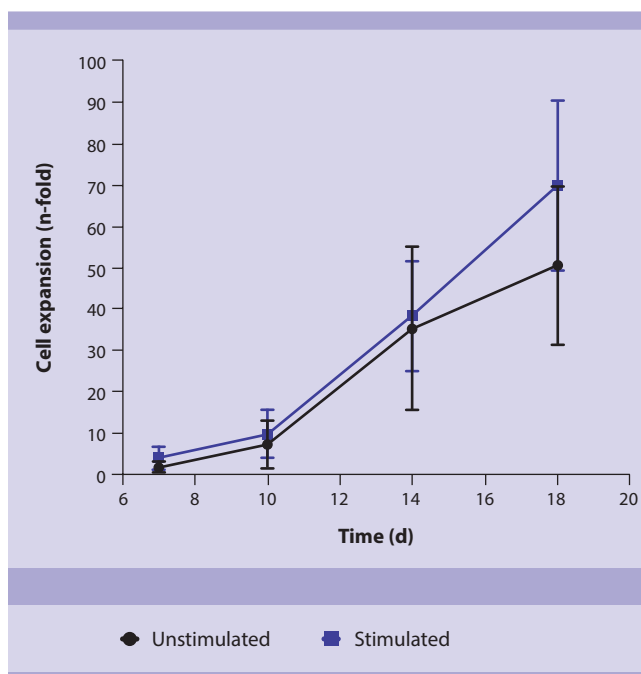

Refer to www.miltenyibiotec.com for all data sheets and protocols. Miltenyi Biotec provides technical support worldwide. Visit www.miltenyibiotec.com for local Miltenyi Biotec Technical Support contact information.

Legal notices

Limited product warranty

Miltenyi Biotec B.V. & Co. KG and/or its affiliate(s) warrant this product to be free from material defects in workmanship and materials and to conform substantially with Miltenyi Biotec's published specifications for the product at the time of order, under normal use and conditions in accordance with its applicable documentation, for a period beginning on the date of delivery of the product by Miltenyi Biotec or its authorized distributor and ending on the expiration date of the product's applicable shelf life stated on the product label, packaging or documentation (as applicable) or, in the absence thereof, ONE (1) YEAR from date of delivery ("Product Warranty"). Miltenyi Biotec's Product Warranty is provided subject to the warranty terms as set forth in Miltenyi Biotec's General Terms and Conditions for the Sale of Products and Services available on Miltenyi Biotec's website at www.miltenyibiotec.com, as in effect at the time of order ("Product Warranty"). Additional terms may apply. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING IF A PRODUCT IS SUITABLE FOR CUSTOMER'S PARTICULAR PURPOSE AND APPLICATION METHODS.

Technical information

The technical information, data, protocols, and other statements provided by Miltenyi Biotec in this document are based on information, tests, or experience which Miltenyi Biotec believes to be reliable, but the accuracy or completeness of such information is not guaranteed. Such technical information and data are intended for persons with knowledge and technical skills sufficient to assess and apply their own informed judgment to the information. Miltenyi Biotec shall not be liable for any technical or editorial errors or omissions contained herein.

All information and specifications are subject to change without prior notice. Please contact Miltenyi Biotec Technical Support or visit www.miltenyibiotec.com for the most up-to-date information on Miltenyi Biotec products.

Licenses

This product and/or its use may be covered by one or more pending or issued patents and/or may have certain limitations. Certain uses may be excluded by separate terms and conditions. Please contact your local Miltenyi Biotec representative or visit Miltenyi Biotec's website at www.miltenyibiotec.com for more information.

The purchase of this product conveys to the customer the non-transferable right to use the purchased amount of the product in research conducted by the customer (whether the customer is an academic or for-profit entity). This product may not be further sold. Additional terms and conditions (including the terms of a Limited Use Label License) may apply.

CUSTOMER'S USE OF THIS PRODUCT MAY REQUIRE ADDITIONAL LICENSES DEPENDING ON THE SPECIFIC APPLICATION. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING FOR ITSELF WHETHER IT HAS ALL APPROPRIATE LICENSES IN PLACE. Miltenyi Biotec provides no warranty that customer's use of this product does not and will not infringe intellectual property rights owned by a third party. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

Trademarks

autoMACS, MACS, MACSiBead, MACSmix, MidiMACS, the Miltenyi Biotec logo, MiniMACS, OctoMACS, QuadroMACS, and SuperMACS are registered trademarks or trademarks of Miltenyi Biotec B.V. & Co. KG and/or its affiliates in various countries worldwide. All other trademarks mentioned in this publication are the property of their respective owners and are used for identification purposes only.

Copyright © 2024 Miltenyi Biotec and/or its affiliates. All rights reserved.