Miltenyi Biotec

gentleMACS™ Octo Dissociator with Heaters

User manual

Copyright © 2026 Miltenyi Biotec and/or its affiliates. All rights reserved.

No part of this user manual may be reproduced, stored in a retrieval system, transmitted, published, or distributed in any form or by any means, electronically, mechanically, by photocopying, microfilming, recording, or otherwise, without the prior written consent of Miltenyi Biotec; however, notwithstanding the foregoing, the owners of the gentleMACS Octo Dissociator with Heaters may make copies solely for purposes of training personnel in the use and servicing of the unit within their business or organization. Maximal care has been taken by Miltenyi Biotec in the preparation of this user manual. However, Miltenyi Biotec shall not be liable for any technical or editorial errors or omissions contained herein, or for incidental or consequential damages in connection with the furnishing, performance, or use of this document. The information in this document is provided "as is" without warranty of any kind, and is subject to change without notice. Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research use only and not for therapeutic or diagnostic use.

gentleMACS™ Octo Dissociator with Heaters

User manual

Original instructions

Issued: 2026-02

140-006-564-006.06

Miltenyi Biotec B.V. & Co. KG
Friedrich-Ebert-Straße 68
51429 Bergisch Gladbach
Germany
Phone +49 2204 8306-0
Fax +49 2204 85197
macsde@miltenyi.com
www.miltenyibiotec.com

Read the chapter **Important safety information** and all other information in this user manual before using the instrument

 WARNING

Before using the gentleMACS Octo Dissociator with Heaters, read the chapter **Important safety information** and all other information contained in this user manual, including any safety and operating instructions. Pay special attention to all warnings displayed on the instrument. Failure to read and follow these guidelines could lead to improper or incorrect usage and result in damage to the instrument. Improper usage could also cause severe personal injury, death, unpredictable results, instrument malfunction, and premature wear to components shortening the lifetime of the instrument. Such actions may void your warranty. Keep the user manual and any other safety and operating instructions provided with the instrument in a safe place accessible to all users for future reference.

If you have a serious concern regarding the safe use of your instrument, contact your authorized Miltenyi Biotec service provider or call Miltenyi Biotec Technical Support.

Supported instruments

This user manual applies to the gentleMACS Octo Dissociator with Heaters (# 130-134-029) running with the gentleMACS Software version 1.2.

Contents

Important safety information	9
Wichtige Sicherheitsinformationen	17
Información importante de seguridad	25
Informations de sécurité importantes	33
Importanti istruzioni di sicurezza	41
1 Introduction	49
1.1 Intended use	49
1.2 Instrument schematics	49
1.3 Sample preparation	50
1.4 gentleMACS Tubes	51
1.5 gentleMACS Heaters	52
1.6 Perfusion Technology	53
1.7 gentleMACS Octo Coolers	53
1.8 MACS Tissue Dissociation Kits	53
1.9 Predefined tissue dissociation and homogenization programs	53
1.10 Customized dissociation programs	53
2 Installation	55
2.1 Components included in the delivery	55
2.2 Requirements	55
2.3 Installation	56
3 Required materials	57
3.1 Disposables	57
3.1.1 Additional disposables	58
3.2 Accessories	58
3.3 Spare parts	58
4 Instrument operation	59
4.1 Switching on	59

4.2 Switching off	59
4.3 The user interface	60
4.4 Home hub	61
4.4.1 Status of position boxes	62
4.5 Programs hub	63
4.5.1 Creating a new program	63
4.5.2 Editing a program	64
4.5.3 Filtering programs	64
4.5.4 Deselecting program filters	65
4.5.5 Deleting programs	65
4.5.6 Exporting programs	65
4.5.7 Importing programs	65
4.6 Settings hub	66
4.6.1 Editing instrument name	66
4.6.2 Editing date and time	67
4.6.3 User settings	67
4.6.4 Instrument settings	68
4.6.5 Software update	68
4.7 LEDs	69
5 Performing runs	71
5.1 Running a single sample	71
5.2 Running multiple samples simultaneously	71
5.3 Running a perfusion program	72
5.4 Pausing and resuming samples	72
5.5 Aborting a running sample	73
6 Maintenance	75
6.1 Decontamination and cleaning	75
7 Troubleshooting	77
7.1 Troubleshooting table	77

7.2 Exporting log files	78
8 Technical support	79
9 Technical data and specifications	81
9.1 FCC compliance statement	82
9.2 EC/EU Declaration of Conformity	83
9.3 UK Declaration of Conformity	84
10 Legal notes	85
10.1 Limited warranty	85
10.2 Trademarks	86
11 Index	87

General safety instructions

THIS CHAPTER DESCRIBES POTENTIALLY HAZARDOUS SITUATIONS ASSOCIATED WITH THIS INSTRUMENT AND PROVIDES IMPORTANT SAFETY INFORMATION TO MINIMIZE THE RISKS AND PROTECT YOURSELF AND OTHERS.

This chapter provides important information for your personal safety and the correct use of the instrument. Read and observe all instructions carefully before proceeding with the installation and use of the instrument. Observe general safety practices in addition to this user manual.

- Use this instrument only as indicated in this user manual to avoid personal injury and property damage.
- Keep this user manual in a place that is always accessible to all users.
- Observe local working area safety instructions and laboratory policies, as well as standards for health, safety, and prevention of accidents.
- In case of severe accidents, damages to the instrument, or if smoke or flames appear, cut the power supply immediately.
- To entirely disconnect the instrument from the power supply, unplug the power cable.
- Ensure that the power switch and the plug of the power cable of the instrument are easily accessible.

Hazard levels

Signal words are used to warn against hazardous situations and property damages. The following signal words are used in this user manual.

⚠ WARNING or **WARNING!** indicates a hazardous situation that, if not avoided, could result in death or serious injury.

⚠ CAUTION or **CAUTION!** indicates a hazardous situation that, if not avoided, could result in minor or moderate injury. It is also used to warn against unsafe practices.

NOTICE or **NOTICE** indicates information considered important but not hazard related (e.g. messages relating to property damage).

Symbols

SAFETY WARNING: THE DOCUMENTATION MUST BE CONSULTED IN ALL CASES WHERE THIS SAFETY SYMBOL IS USED ON THE INSTRUMENT

HOT SURFACE

ROTATING PARTS

ON (POWER ON)

OFF (POWER OFF)

READ THE USER MANUAL BEFORE USING THE INSTRUMENT

FUSE

WEEE (WASTE ELECTRICAL AND ELECTRONIC EQUIPMENT)

ORDER NUMBER

TYPE NUMBER

SERIAL NUMBER

MANUFACTURER

DATE OF MANUFACTURE

EUROPEAN CONFORMITY MARKING


UNITED KINGDOM CONFORMITY ASSESSED MARKING

NRTL CERTIFICATION MARK: PRODUCT MEETS CONSENSUS-BASED STANDARDS OF SAFETY, REQUIRED BY THE OCCUPATIONAL SAFETY/HEALTH ADMINISTRATION (OSHA), DETERMINED BY THE NATIONALLY RECOGNIZED TESTING LABORATORIES (NRTL) TÜV SÜD

Safety labels

Notice the hazard points and safety symbols of the instrument. Keep safety labels and safety markings clean and legible. Inspect the safety labels and safety markings regularly and replace them if they are not legible or identifiable from a safe viewing distance. Contact Miltenyi Biotec for replacement labels. The following safety label is attached to the back of the instrument and plays an important part in maintaining safety.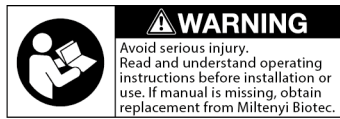

WARNING! Avoid serious injury. Read and understand operating instructions before installing or use. If manual is missing, obtain replacement from Miltenyi Biotec.

The gentleMACS Sleeves are marked with a contrasting safety print. The contrasting safety print allows the visual perception of the rotation of the gentleMACS Sleeves. The gentleMACS Octo Trough is marked with the following safety symbol.

Figure 1: Safety symbol on the gentleMACS Octo Trough and contrasting safety print on the gentleMACS Sleeves

- 1 Contrasting safety print on the gentleMACS Sleeves
- 2 Safety symbol on the gentleMACS Octo Trough

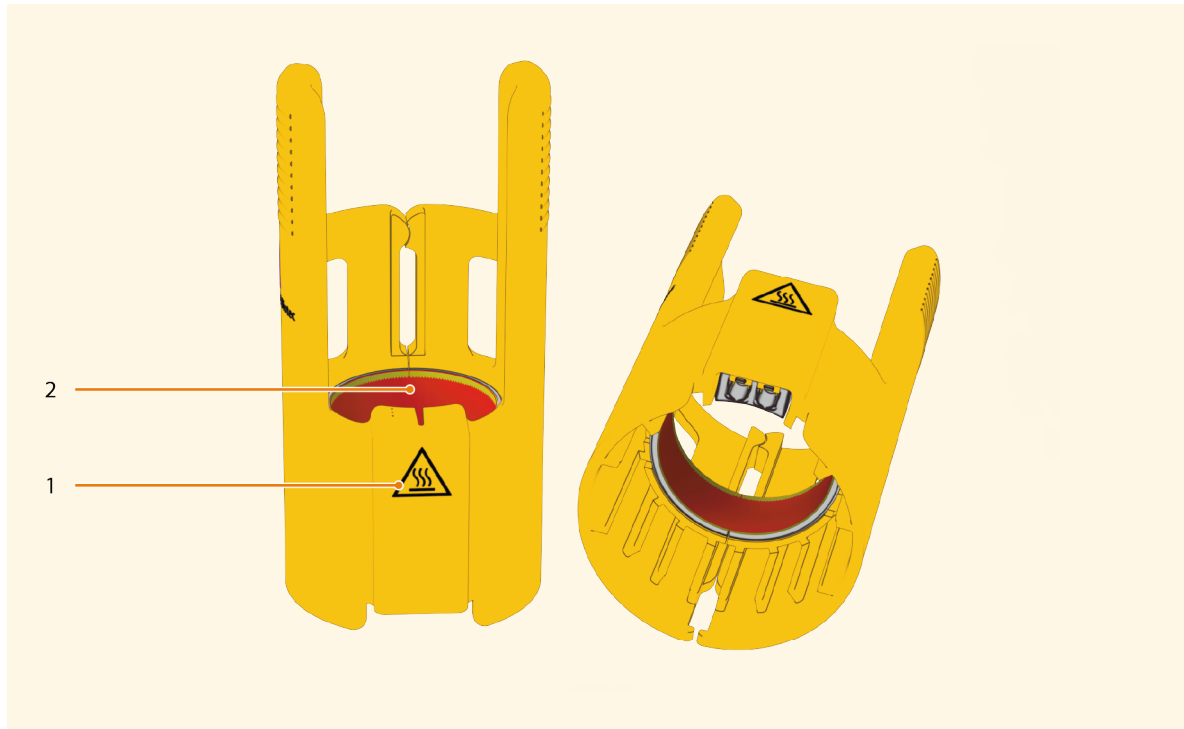
Figure 2: Safety symbol on the gentleMACS Heaters

- 1 Safety symbol on the gentleMACS Heaters
- 2 Heating foil

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Protective housing

The housing of the instrument reduces the risk of an electric shock and short circuits.

- Do not remove or penetrate any cover of the housing.
- Exceptions are maintenance procedures that are explicitly described in this manual.
- Do not use the instrument if it is opened or disassembled.
- Do not use the instrument if an object entered the inside of the instrument.
- Check the instrument before each use for visible damages.
- Do not use the instrument if it has been dropped or is damaged.

Cables and power supplies

Using the instrument with other than the supplied cables is potentially hazardous.

- Only use the included power cable.
- Do not use the instrument if the power cable is damaged.
- The instrument has a three-wire electrical grounding plug with a third pin for grounding.
- Only insert the plug into a grounded electrical outlet. Do not try to insert the plug into a non-grounded electrical outlet.
- Operate the instrument only from a power source that meets the electrical specifications mentioned on the type label. If you have questions about the type of power source to use, contact the Miltenyi Biotec Technical Support or your local electricity supplier.

Liquids inside the instrument

Liquids inside the instrument can cause short circuits.

- Protect the instrument against spilled liquids.
- Do not use the instrument if liquids have entered the inside of the instrument.
- If liquids have entered the instrument, unplug the instrument and contact Miltenyi Biotec Technical Support.
- Clean up spilled liquids immediately.
- Unplug the instrument before cleaning it.
- Use only small amounts of cleaning agents on a soft cloth to wipe the instrument. Do not spray or pour liquid cleaning agents onto or into the instrument.

Mechanical hazards

CAUTION

Risk of crushing.

The instrument will fall down if it is not placed on a flat and even surface, which could result in injuries.

- Place the instrument only onto stable, even, and vibration free tables or laboratory benches.
- Handle the instrument with care during carrying or relocation.

CAUTION

Risk of crushing or cutting.

gentleMACS Tube caps have sharp edges on the inside. This may lead to incision wounds if they were touched while rotating.

- Only use tightly closed tubes and ensure that they are properly installed.
- Do not install or remove tubes while the rotors of the instrument are in motion.

CAUTION

Risk of trapping fingers and hair pulled in.

Hair can be pulled in by the rotors when the instrument is running.

- Do not operate the instrument without gentleMACS Octo Trough, gentleMACS Sleeves, and gentleMACS Octo Safety Shield.
- Tie back long hair.
- Do not touch the rotors while the instrument is running.

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Biological material

All biological material must be considered potentially infectious.

- Wear personal protective equipment (such as gloves, safety glasses, etc.) as indicated in the safety data sheet for the particular substance.
- Operate the instrument in a safety hood if hazardous or potentially infectious materials are processed.
- Decontaminate the instrument after processing hazardous or potentially infectious material.
- If hazardous or potentially infectious material has been spilled or leaked from the system, decontaminate the affected area.
- Use a disinfectant or cleaning agent that is appropriate for the specific pathogen or substance.

Waste

All liquid and solid waste must be considered hazardous.

- Observe local regulations regarding waste disposal.
- Autoclave or alternatively decontaminate waste with an appropriate disinfectant.

Chemical hazards

WARNING

Substances and reagents can be hazardous.

- All safety measures in section **Biological hazards** also apply to any hazardous substances and reagents that may be present in the sample.
- Operate the instrument in a fume hood if hazardous substances and reagents are processed.

Thermal hazards

CAUTION

Hot surfaces can cause burns.

When the heating function is used, the heating foil of the gentleMACS Heaters becomes hot.

- Do not touch the heating foil.

Noise

CAUTION

High noise levels could result in hearing impairment.

If running with maximum load, the instrument may produce loud noise. It is recommended that the responsible body measures or calculates the sound pressure level at the operator's position and at whatever point 1 m from the enclosure of the equipment has the highest sound pressure level.

- Noise level should be below 70 dB.
- Observe local regulations regarding hearing protection.
- If running the instrument with maximum load, then protective safety equipment such as noise reducing baffles or hoods, or ear protection has to be used.

Servicing and transportation

Servicing

Improper servicing or repair of the instrument or use of unauthorized parts can cause malfunction of or damage to the instrument. This can cause hazards to users.

- Unless otherwise specifically stated in this user manual or other documentation by Miltenyi Biotec, do not service the instrument yourself.
- Servicing and repair must be performed by Miltenyi Biotec certified and qualified service personnel.
- If the instrument needs servicing, decontaminate the instrument to remove any hazardous material.
- If you have questions regarding proper decontamination or shipment, contact Miltenyi Biotec Technical Support for assistance.
- Only use accessories and upgrades recommended by Miltenyi Biotec.
- Inquire with your local Miltenyi Biotec representative about Miltenyi Biotec's extensive instrument service and service contracts, or refer to **www.miltenyibiotec.com/support**.

Transportation

The instrument must be transported with care in packaging specified by Miltenyi Biotec. Internal damage could occur if the instrument is subjected to excessive vibration or if it is dropped. If the instrument needs to be shipped back to the manufacturer for service, contact Miltenyi Biotec for instructions and packaging materials.

Disposal

WASTE ELECTRICAL AND ELECTRONIC EQUIPMENT (WEEE) CUSTOMER INFORMATION

Dispose of your end-of-life Miltenyi Biotec products in accordance with the applicable WEEE and hazardous waste disposal legislation, which may differ by country or region. Electrical equipment may contain hazardous substances that may have a serious detrimental effect on the environment and/or human health. All equipment must be specifically collected and treated by designated waste facilities and by qualified WEEE compliance schemes. By ensuring that you dispose of your unwanted electrical and electronic equipment according to the applicable WEEE and hazardous waste disposal legislation, you are helping to preserve our natural resources and protect human health. Miltenyi Biotec is committed to protecting the environment. Miltenyi Biotec offers product end-of-life return programs in many countries and partners with licensed WEEE compliance schemes throughout the world. Miltenyi Biotec takes back your end-of-life Miltenyi Biotec equipment for recycling free of charge. The terms and availability of this offer vary by geography because of differences in regulatory requirements. Note that, depending on the type and use of your equipment, additional requirements may apply. Before shipping the instrument back to the manufacturer for disposal, decontaminate the instrument to remove any hazardous material. For more information or if you wish to dispose of your end-of-life Miltenyi Biotec equipment, contact your local Miltenyi Biotec representative or Miltenyi Biotec Technical Support.

EN				
DE				
ES				
FR				
IT				

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1.1 Intended use

The gentleMACS Octo Dissociator with Heaters is a benchtop instrument for fully automated and standardized tissue dissociation, perfusion or homogenization of up to eight samples. Equipped with eight individual heating units and numerous optimized and ready-to-use gentleMACS Programs, the gentleMACS Octo Dissociator with Heaters offers a fully automated workflow for tissue dissociation, perfusion, and homogenization. The gentleMACS Octo Dissociator with Heaters is for research-use only in a laboratory environment. It is to be used by laboratory professional users only.

1.2 Instrument schematics

- | | | | |
|---|--|----|---|
| 1 | gentleMACS Safety Shield, referred to as Safety Shield | 7 | gentleMACS Sleeve, referred to as Sleeve |
| 2 | Touchscreen | 8 | LAN port |
| 3 | Power switch | 9 | CAN port |
| 4 | Power port | 10 | Type label |
| 5 | USB port | 11 | gentleMACS Octo Trough, referred to as Trough |
| 6 | 2D barcode reader | | |

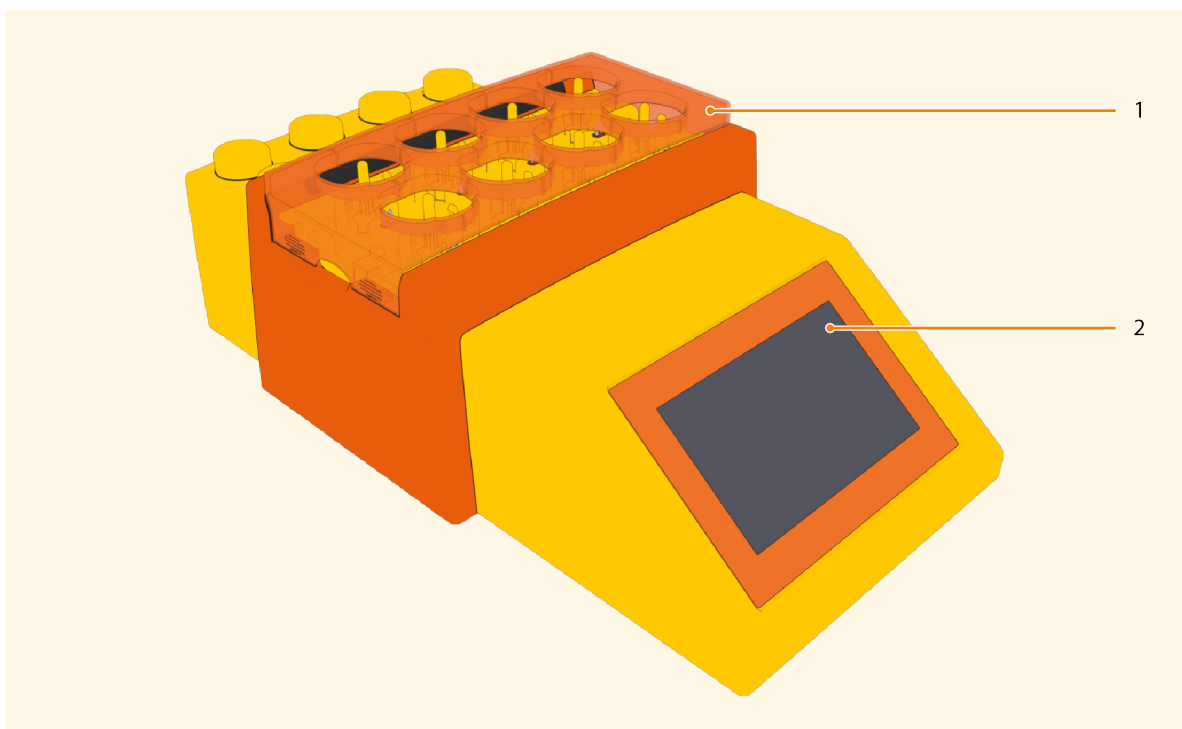
Figure 1.1: Front view of the gentleMACS Octo Dissociator with Heaters

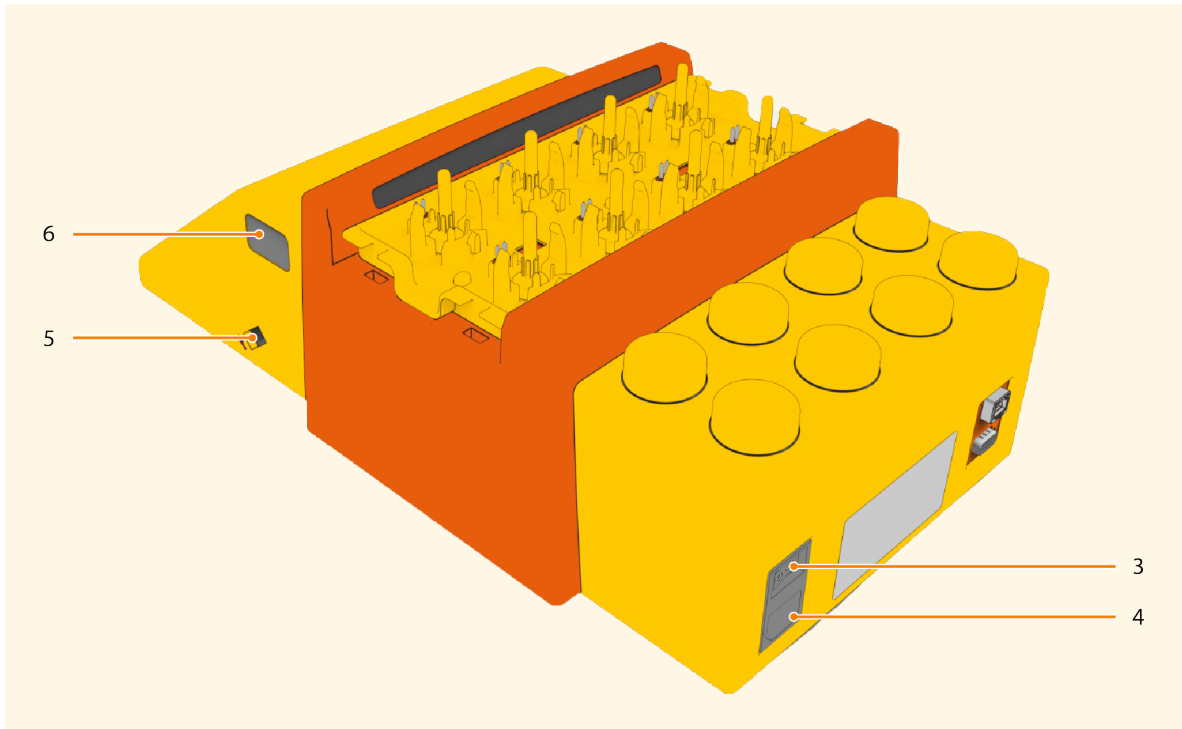
Figure 1.2: Rear view of the gentleMACS Octo Dissociator with Heaters, Safety Shield removed

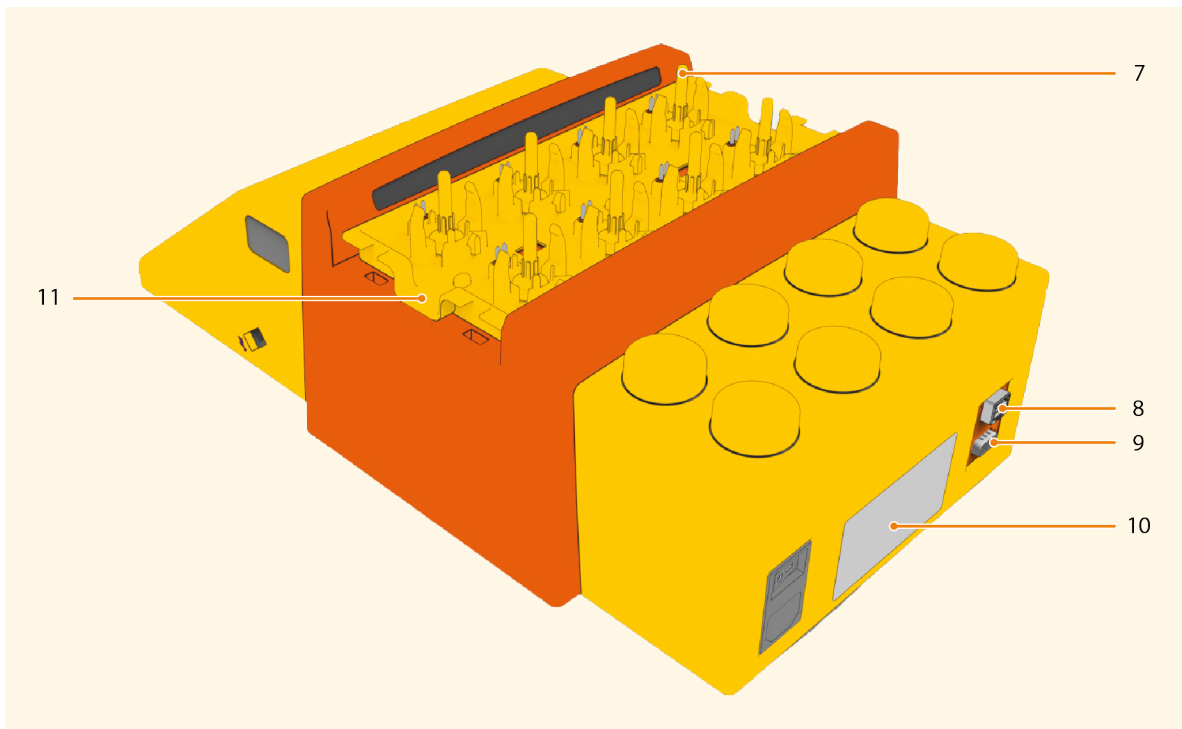
Figure 1.3: Rear view of the gentleMACS Octo Dissociator with Heaters, Safety Shield removed

1.3 Sample preparation

The gentleMACS Octo Dissociator with Heaters provides the solution for the preparation of single-cell, or single-nuclei suspensions, and homogenates from a wide range of tissues. Samples are processed using gentleMACS Tubes, which have caps with specially designed rotors and stators.

The gentleMACS Octo Dissociator with Heaters provides:

- the preparation of viable single-cell suspensions by mechanical and enzymatic dissociation, including incubations at 37 °C and in combination with the gentleMACS Octo Coolers also between +2°C to +8°C,
- the preparation of intact nuclei suspension by mechanical disruption at low temperature in combination with the gentleMACS Octo Coolers (+2°C to +8 °C),
- the preparation of thorough tissue homogenates for subsequent extraction of sub-cellular material, such as total RNA, DNA, proteins, organelles, and pathogens,
- the *ex vivo* perfusion of resected rodent organs, such as liver and heart, using the gentleMACS Perfusion Technology,
- simultaneous or independent processing of up to eight samples, even using different programs,
- easy access to predefined gentleMACS Programs which are optimized for particular applications,
- the option to create user-defined programs.

1.4 gentleMACS Tubes

gentleMACS Tubes consist of a stator and rotor element that provides the exact sheer forces necessary to gently pull cells from tissues or intact molecules from cells. A specially engineered tube enclosure always directs the sample flow towards the rotor and stator element. This sophisticated design results in highly efficient tissue dissociation and homogenization using C Tubes and M Tubes.

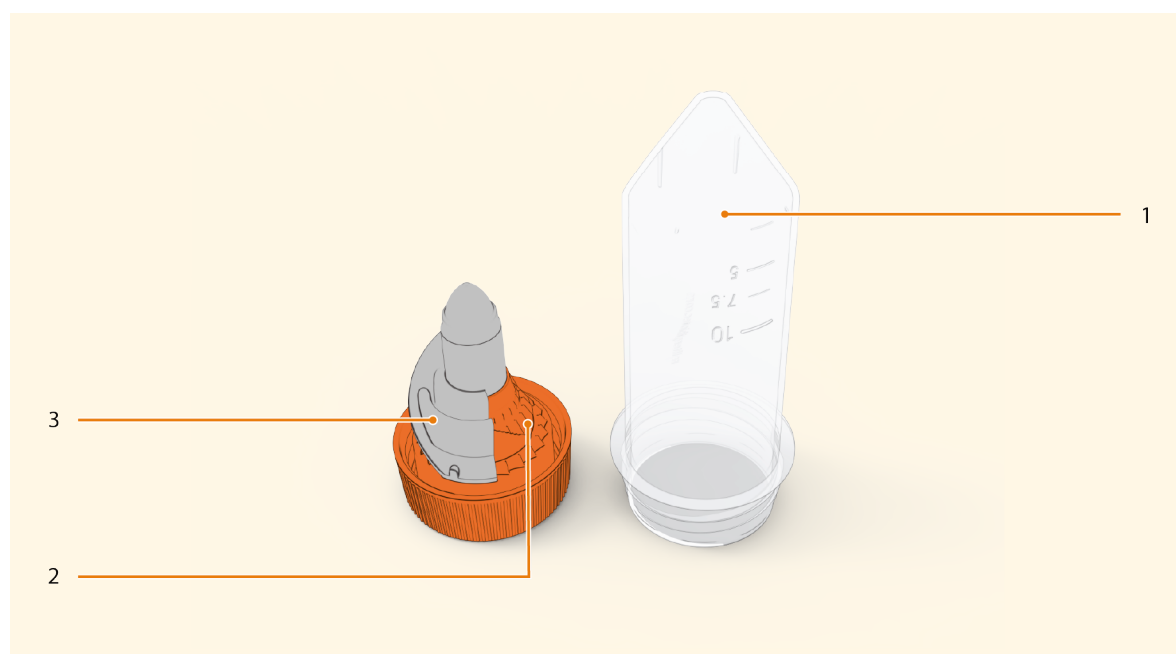
Figure 1.4: Cap of the gentleMACS Tubes (exemplary)

1 Tube enclosure

3 Rotor

2 Stator

Two dedicated types of gentleMACS Tubes have been designed for the preparation of samples for cell biology or molecular biology applications:

- Purple C Tubes are designed for the gentle preparation of single-cell, or single-nuclei suspensions from tissues for subsequent cell separation, cell analysis, cell culture, or genomic analysis, respectively.
- Orange M Tubes are designed for the homogenization of tissues and cells for subsequent subcellular or molecular analysis.

gentleMACS Tubes are suitable for:

- the processing of sample volumes from 500 μ L to 10 mL and from 20 mg to 4,000 mg of tissue in a single tube,
- sterile sample handling.

1.5 gentleMACS Heaters

For combined enzymatic and mechanical dissociation processes the instrument is equipped with gentleMACS Heaters for each of the eight tube positions. The gentleMACS Heaters allow to apply 37 °C either to all eight positions or individually to each tube. By gently pressing the clamps, the gentleMACS Heaters can be easily attached to the tube positions, thus connecting the gentleMACS Heaters to the instrument via its electrical contact pins.

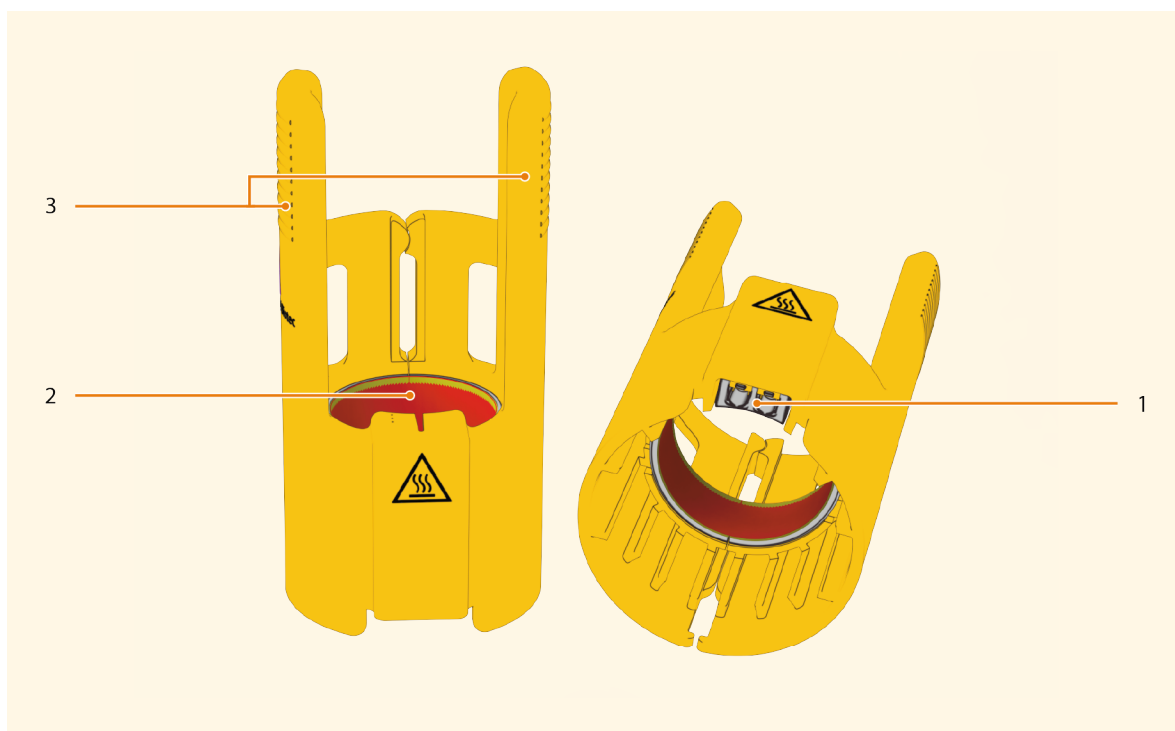
Figure 1.5: The gentleMACS Heaters for incubation of samples at 37 °C

1 Electrical contacts

3 Clamps

2 Heating foil

1.6 Perfusion Technology

The gentleMACS Perfusers are designed for the gentleMACS Perfusion Technology and are used in combination with the gentleMACS Perfusion Sleeves and the gentleMACS Octo Dissociator with Heaters. They are single-use consumables that enable *ex vivo* perfusion of tissues. During the perfusion process, enzymes digest the tissue. Subsequent processing of the tissue with C Tubes results in a single-cell suspension which can be used for a wide variety of different downstream applications, such as cell separation, cell culture, cell or molecular analysis.

1.7 gentleMACS Octo Coolers

The gentleMACS Octo Coolers are an accessory for gentleMACS Octo Dissociators. They are designed to maintain the temperature of samples inside the tubes between +2°C to +8 °C during cold dissociation or homogenization. The coolers are compatible with all types of gentleMACS Tubes or gentleMACS Perfusers.

1.8 MACS Tissue Dissociation Kits

A broad selection of MACS Tissue Dissociation Kits, including predefined enzyme mixes, provide convenient and reproducible tissue dissociation using the gentleMACS Octo Dissociator with Heaters. A comprehensive list of MACS Tissue Dissociation Kits for numerous applications is available for download at www.miltenyi.com/epitopes.

1.9 Predefined tissue dissociation and homogenization programs

More than 80 preset gentleMACS Programs have been specifically developed to offer standardized procedures for tissue dissociation in virtually any research field. A comprehensive list of optimized gentleMACS Programs and detailed gentleMACS Protocols for numerous applications are available for download at www.miltenyi.com/brochureapps.

1.10 Customized dissociation programs

In addition to the ready-to-use gentleMACS Programs optimized for specific applications, the gentleMACS Octo Dissociator with Heaters also allows users to create new programs. See **Creating a new program on page 63**.

2

Installation

Unpack the instrument carefully and confirm that all components have been successfully delivered without damage that may have occurred during transportation.

WARNING! If there is any damage, do not use the instrument. Contact your local Miltenyi Biotec representative or Miltenyi Biotec Technical Support. Operating a damaged instrument may cause mechanical hazards, optical radiation hazards, fire hazards, or electric shock. Read the chapter Important safety information before installing and assembling the instrument.

Retain the original packaging material.

2.1 Components included in the delivery

The following components are included in the delivery of the gentleMACS Octo Dissociator with Heaters (# 130-134-029):

- 1× gentleMACS Octo Dissociator with Heaters
- 8× gentleMACS Heaters
- 1× Power cable
- 1× User manual

2.2 Requirements

Meet the following requirements to install the instrument:

- The instrument requires a power supply.
- Place the instrument on a flat and clean surface of the dimensions 495 mm × 285 mm plus at least 15 cm on all sides to ensure for sufficient air circulation.
- The display and top of the instrument must be accessible.
- The instrument must be operated in a safety hood if hazardous or unknown materials are being processed.
- Do not place the instrument near radiators, heat registers, stoves, or other equipment (including amplifiers) that generate heat.
- Protect the instrument from direct sunlight.

2.3 Installation

To install the gentleMACS Octo Dissociator with Heaters:

- 1 Carefully remove the device from the cardboard box.

CAUTION! The instrument could fall down and cause serious injury to hand or foot. Place the instrument only onto stable, even, and vibration free tables or laboratory benches.

- 2 Place the instrument on a secure, clean, and stable workbench that can support its weight and size.
- 3 Plug the instrument into an electrical outlet.
- 4 Set the switch located above the plug socket to ON ("I" indicates "ON", "O" indicates "OFF").

The instrument performs a start-up check that runs for a few seconds.

3

Required materials

3.1 Disposables

The materials listed in **Table 3.1** are compatible with the gentleMACS Octo Dissociator with Heaters.

Product	Description	Capacity	Order no.
gentleMACS C Tubes	For gentle preparation of - single-cell suspensions from tissues for subsequent cell separation, cell analysis, and cell culture. - intact nuclei suspension for genomic analysis.	25 tubes, sterile, single-packed	130-093-237
		100 tubes, sterile, packed as 4x25 tubes	130-096-334
gentleMACS M Tubes	For homogenization of tissues and cells for subsequent subcellular and molecular analysis.	25 tubes, sterile, single-packed	130-093-236
		100 tubes, sterile, packed as 4x25 tubes	130-096-335
M Tubes with Strainer	For homogenization of tissues and cells for subsequent subcellular and molecular analysis. The strainer removes aggregates or large particles from the sample.	50 tubes, sterile, packed as 2x25 tubes (strainer: 600 µm mesh)	130-094-392
gentleMACS Perfusers	For <i>ex vivo</i> perfusion of tissues, in combination with the gentleMACS Perfusion Sleeves.	25 tubes, sterile, single-packed	130-128-151
gentleMACS Perfusers 2	For <i>ex vivo</i> perfusion of smaller tissues, in combination with the gentleMACS Perfusion Sleeves.	25 tubes, sterile, single-packed	130-134-803

Table 3.1: Disposables compatible with the gentleMACS Octo Dissociator with Heaters

3.1.1 Additional disposables

MACS SmartStrainers are filters that are designed for the removal of cell aggregates or large particles after tissue dissociation. The MACS SmartStrainers are available with three mesh sizes:

- 30 µm (50 pieces # 130-098-458, 4×25 pieces # 130-110-915),
- 70 µm (50 pieces # 130-098-462, 4×25 pieces # 130-110-916),
- 100 µm (50 pieces # 130-098-463, 4×25 pieces # 130-110-917).

They fit into standard 15 and 50 mL conical tubes of all major suppliers.

MACS Tissue Dissociation Kits were designed for the convenient and standardized enzymatic dissociation of multiple tissues in combination with the gentleMACS Octo Dissociator with Heaters. For a comprehensive list of MACS Tissue Dissociation Kits, see www.miltenyi.com/epitopes.

3.2 Accessories

Product	Description	Order no.
gentleMACS Octo Coolers (4×)	For protocols with a cooling step	130-130-533
gentleMACS Perfusion Sleeves (sterile, packed as 2×2 pieces)	Adapters for attaching gentleMACS Perfusers and gentleMACS Perfusers 2	130-128-752

Table 3.2: Accessories for the gentleMACS Octo Dissociator with Heaters

3.3 Spare parts

Product	Description	Order no.
gentleMACS Sleeves (2×)	Adapters for attaching gentleMACS Tubes	130-093-464
gentleMACS Heater	Heating unit	130-105-228
gentleMACS Octo Cooler (1×)	For protocols with a cooling step	130-131-248
gentleMACS Octo Trough	Safety container for spilled liquid	130-136-426
gentleMACS Octo Safety Shield	Safety shield for protection	130-096-542

Table 3.3: Spare parts of the gentleMACS Octo Dissociator with Heaters

Only use spare parts and accessories recommended by Miltenyi Biotec. Failure to use recommended spare parts and disposables may compromise safe operation and may result in inaccurate results and reduced lifetime of the instrument. Miltenyi Biotec does not honor any warranty or accept any responsibility for damages resulting from the use of inappropriate accessories.

4

Instrument operation

4.1 Switching on

Switch on the instrument with the power switch on the rear side of the instrument (**Figure 1.2**).

The instrument starts and the software opens. The internal fan automatically starts operating and reduces its speed after a few seconds.

After the first start of the instrument, naming the instrument and setting date and time is required. This information can be edited. See **Editing instrument name on page 66** and **Editing date and time on page 67**.

4.2 Switching off

- 1 Shut down the software before switching off the instrument. Tap the shut down button in the top indicator bar (**Figure 4.1**).
- 2 Tap **Shutdown** in the message window to proceed.
The software shuts down.
- 3 Switch off the instrument with the power switch on the rear side of the instrument (**Figure 1.2**).

4.3 The user interface

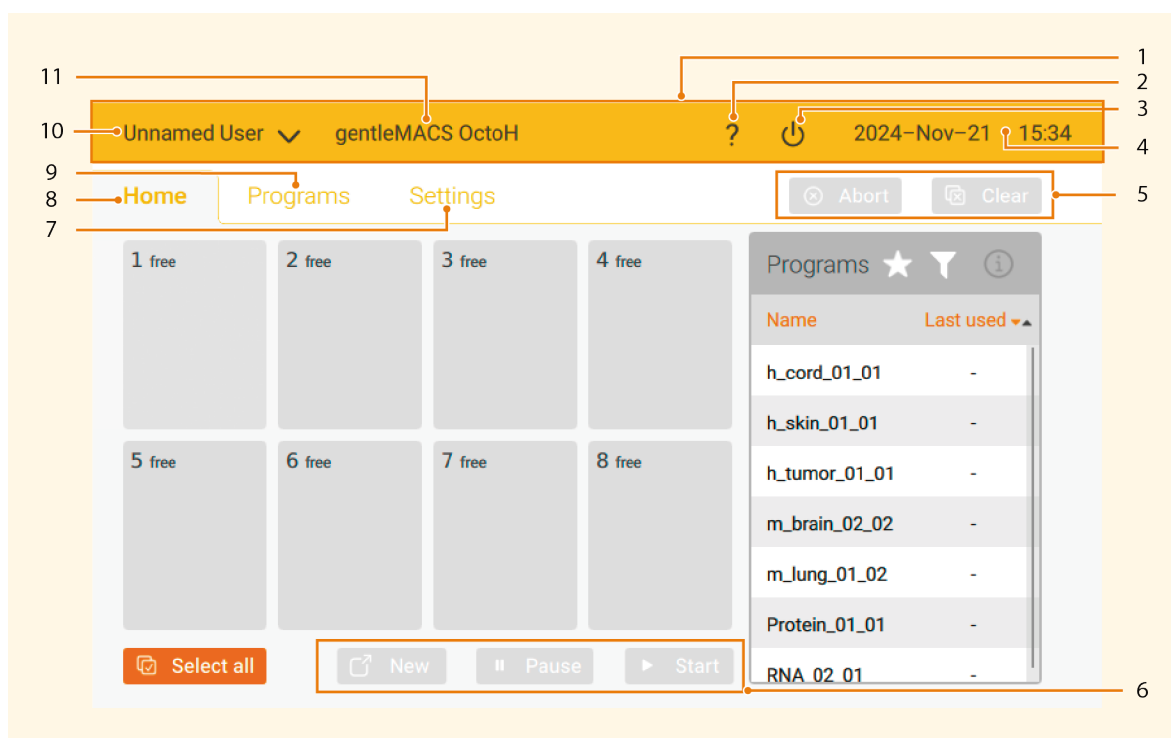
Figure 4.1: Overview of the user interface

1	Top indicator bar	Shows name of current user, instrument name, date and time; includes the help button and shut down button
2	Help button	Opens the help information on the instrument
3	Shut down button	Button to shut down the software before switching off the instrument
4	Date and time	Current date and time set on the instrument
5	Local position buttons	Buttons controlling the functions Abort , and Clear for all selected positions
6	Global position buttons	Buttons controlling the functions New , Pause and Start for all positions in the corresponding status
7	Home hub	Menu hub with instrument position overview and Programs list; gives an overview of the status of each position and is the starting point for assigning programs to positions and controlling sample runs
8	Programs hub	Menu hub for managing programs, including the program editor
9	Settings hub	Menu hub for managing instrument and user settings
10	User	Shows the current user of the instrument, users can be switched in the dropdown list
11	Instrument name	Name of the instrument

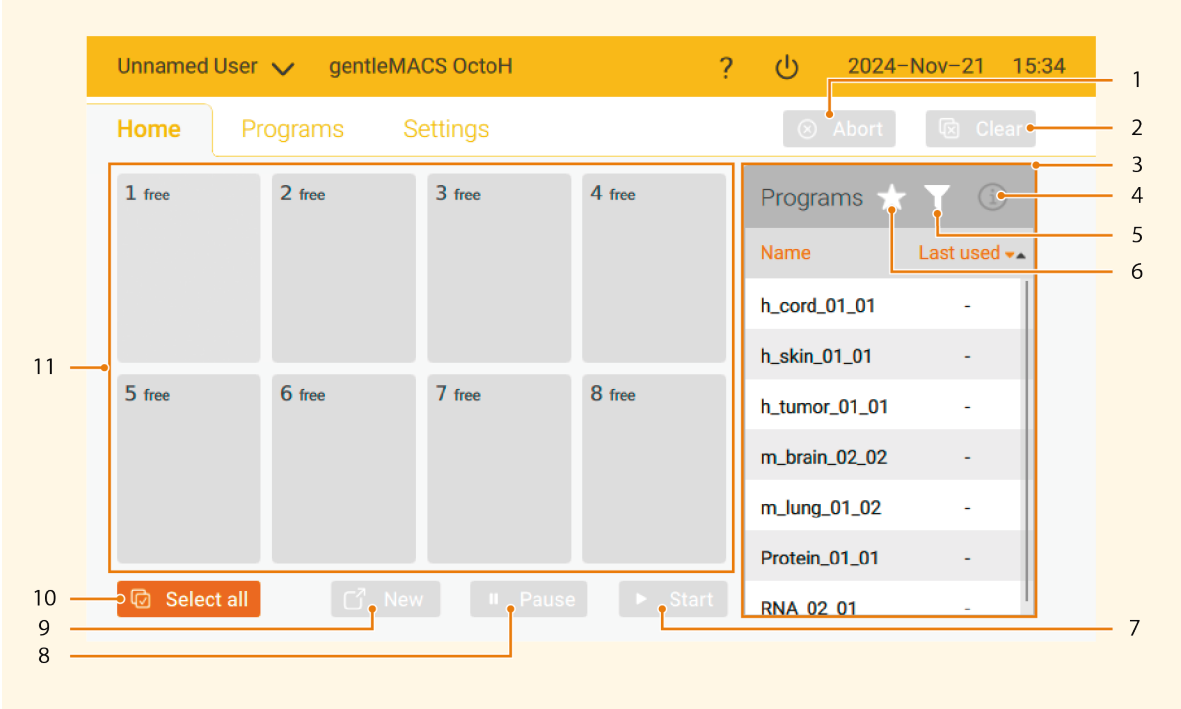
Figure 4.2: Overview Home hub

1	Abort button	Aborts run of the selected position boxes
2	Clear button	Clears the assigned programs or Done status of the selected position boxes
3	Programs list	List of programs available
4	Program details button	Shows information about the selected program like tube type needed, duration, rotations per run, species, starting material etc.; option to mark selected program as Favorite
5	Filter button	Opens Filter attribute window; option to filter by tube type, thermal condition, species, starting material, program author and favorite
6	Favorites button	Filters for favorite programs of the logged-in user
7	Start button	Starts runs on all positions in the status Ready
8	Pause button	Pauses all active sample runs
9	New button	Clears the assigned programs of all position boxes that are not running or paused
10	Select all/ Deselect all button	Selects or deselect all position boxes at once. If all position boxes are selected, the Select all button turns into a Deselect all button.
11	Position boxes	Managing the runs on the eight instrument positions; shows status of each position (see Status of position boxes on the next page)

4.4.1 Status of position boxes

Status	Color	Description
free	Gray	No tube inserted, no program assigned. The position is free to use.
selected	Gray with orange border	The position is selected.
partly	Green, pulsing	Position is partly ready. A program is assigned to the position box, but a tube and/or heater is still needed, indicated by the icon in the upper right corner of the position box. Or a tube and/or heater are placed but no program is assigned to the position box.
ready	Green	The position is ready. A tube and/or heater is inserted and a program is assigned.
running	Blue	A program is running on the position. The progress bar and timer display the remaining time to complete the program.
paused	Blue, pulsing	The current running program on the position is paused.
break	Blue, pulsing	The current running program on the position has reached a break state. To resume the program, tap the local Start button in the position box and tap Resume in the pop-up window.
done	Gray	The program is finished. The position details button shows the elapsed time since completion of the program. The position box status changes to Done also after a run is aborted.
error	Red	An error occurred. See Troubleshooting table on page 77 .

Table 4.1: Overview of position box statuses

4.5 Programs hub

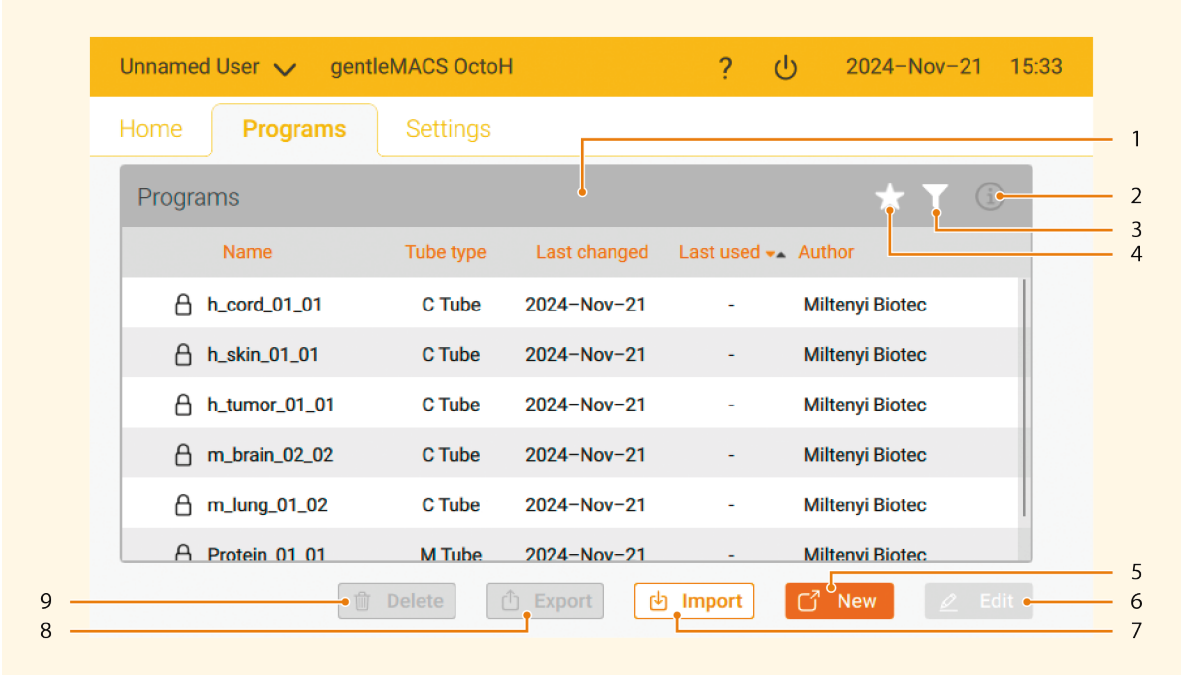
Figure 4.3: Overview Programs hub

- | | | |
|---|------------------------|---|
| 1 | Programs list | List of programs available. The padlock icon indicates which programs can be edited or deleted.
Programs can be sorted by Name, Tube type, Last changed, Last used, and Author by tapping the row heading. |
| 2 | Program details button | Shows information about the selected program like tube type needed, duration, total rotations per run, species, starting material etc.;
option to mark selected program as Favorite . |
| 3 | Filter button | Opens filter attribute window with the option to filter by tube type, thermal condition, species, starting material, program author, and favorite. |
| 4 | Favorites button | Filters for favorite programs of the current user. |
| 5 | New button | Create new program. Opens the program editor window. |
| 6 | Edit button | Edit selected program. Opens the program editor window. Only user programs can be edited. |
| 7 | Import button | Imports selected program from the USB flash drive. |
| 8 | Export button | Exports selected program to USB flash drive. |
| 9 | Delete button | Delete selected program. Only custom programs can be deleted. |

4.5.1 Creating a new program

Do not exceed 3,200 rotations per run, otherwise the gentleMACS Tubes might get damaged.

- 1 Tap **New** in the **Programs** hub. The program editor window opens.
- 2 Enter a name for the new program.
- 3 Tap on the plus symbol to insert a new command. The position where a further command is inserted can be changed using the arrow buttons.
- 4 Add and define all commands needed.
- 5 If needed, commands can be changed, deleted, or rearranged.
Change the position to insert a further command using the arrow buttons if needed.

- 6 Tap **Edit filter** to add filter parameters for the program. At least the tube type needed to run the program has to be defined.
- 7 Tap **Save** to save the new program.
- 8 To return to the **Programs** list, tap the back button next to the program name.

Program commands

Command	Description
Spin	Defines the speed, direction, and duration of rotation in seconds. Speed is defined as revolutions per minute (rpm). The change of velocity must not exceed 6,000 rpm per second.
Heater	Defines if heaters are turned on or off. Add a command to turn heaters on and another to turn heaters off. The temperature of the heaters is set to 37 °C and can not be changed.
Loop	Adds start and end of a loop to the program. Commands between the start and end loop commands are repeatedly executed for the defined numbers of cycles.
Ramp	Defines the time in seconds required to reach a certain speed in rpm. If a program starts with a ramp command, the starting speed will be 0 rpm. If a ramp command is added after a spin command, the starting speed will be from the rpm of the spin command.
No spin	Defines duration in seconds of a step without spinning
Incubate	Defines the duration in minutes of the incubation step with a slow spinning of 20 rpm counterclockwise
Break	Adds a break command. The program will pause in a break state until the user resumes the program.

Table 4.2: Overview of program commands for custom programs

4.5.2 Editing a program

Only custom programs can be edited or deleted. Custom programs do not show the padlock symbol.

- 1 Open the **Programs** hub.
- 2 Tap on the program to be edited.
- 3 Tap **Edit**. The program editor windows opens.
- 4 Edit the program. You can change the program name. Commands can be changed, deleted or rearranged. To change the order of commands, tap on the respective command to select it and change position with the arrow buttons.
- 5 Tap **Edit filter** to change filter parameters for the program. At least the tube type needed to run the program has to be defined.
- 6 Tap **Save** to save the changed program.

4.5.3 Filtering programs

The system combines multiple filter attributes. Attributes of the same category are combined with the filter requirement "or", attributes from different categories are combined with the filter requirement "and".

Users can filter the **Programs** list for programs with certain attributes.

- 1 Tap on the filter button in the **Home** hub or **Programs** hub to open the **Filter** window.

- 2 Select all the attributes that the program should have. Filter attributes are:
 - tube type,
 - thermal condition,
 - species,
 - starting material,
 - author, and
 - favorite.
- 3 Tap X to close the **Filter** window. The **Programs** list shows only programs with the selected attributes.

4.5.4 Deselecting program filters

To remove the filter applied to the program list, perform the following steps:

- 1 Tap on the filter button in the **Home** hub or **Programs** hub to open the **Filter** window.
- 2 Tap **Deselect all**.
- 3 Tap X to close the **Filter** window.

4.5.5 Deleting programs

Only custom programs can be edited or deleted.

Assigned programs can not be deleted. Miltenyi Programs can not be deleted.

- 1 Open the **Programs** hub.
- 2 Tap on the program to be deleted in the programs list.
- 3 Tap **Delete**.
- 4 Confirm the deletion of the selected program.

4.5.6 Exporting programs

You can export programs to a USB flash drive and transfer them to another gentleMACS instrument. Programs are exported as bundles containing one or several programs. Make sure that the USB flash drive is formatted as exFAT for compatibility with the operating system. If required, change the format on a PC.

- 1 Insert the USB flash drive to the USB port of the gentleMACS Octo Dissociator with Heaters.
- 2 Go to the **Programs** hub.
- 3 Tap on the program to export.
To select several programs for export, tap and hold on the first program and then select all other programs to be exported.
- 4 Tap **Export**.
- 5 Choose the target folder on the USB flash drive to export the program to.
- 6 Confirm to resume the export to the USB flash drive.
- 7 After the export is finished, remove the USB flash drive from the instrument.

4.5.7 Importing programs

Use this function to import programs from a USB flash drive to transfer them to another gentleMACS instrument, or to update the Miltenyi Biotec program database with new Miltenyi Biotec programs. Only programs which are supplied by Miltenyi Biotec, with Miltenyi Biotec as specified author, are permitted for

import. Programs are imported as bundles containing one or several programs. Make sure that the USB flash drive is formatted as exFAT for compatibility with the operating system. If required, change the format on a PC.

- 1 Insert the USB flash drive to the USB port of the gentleMACS Octo Dissociator with Heaters.
- 2 Go to the **Programs** hub.
- 3 Tap **Import**.
- 4 Choose the program bundles to be imported from the USB flash drive.
- 5 Confirm to resume the import from the USB flash drive.
- 6 After the import is finished, remove the USB flash drive from the instrument.

4.6 Settings hub

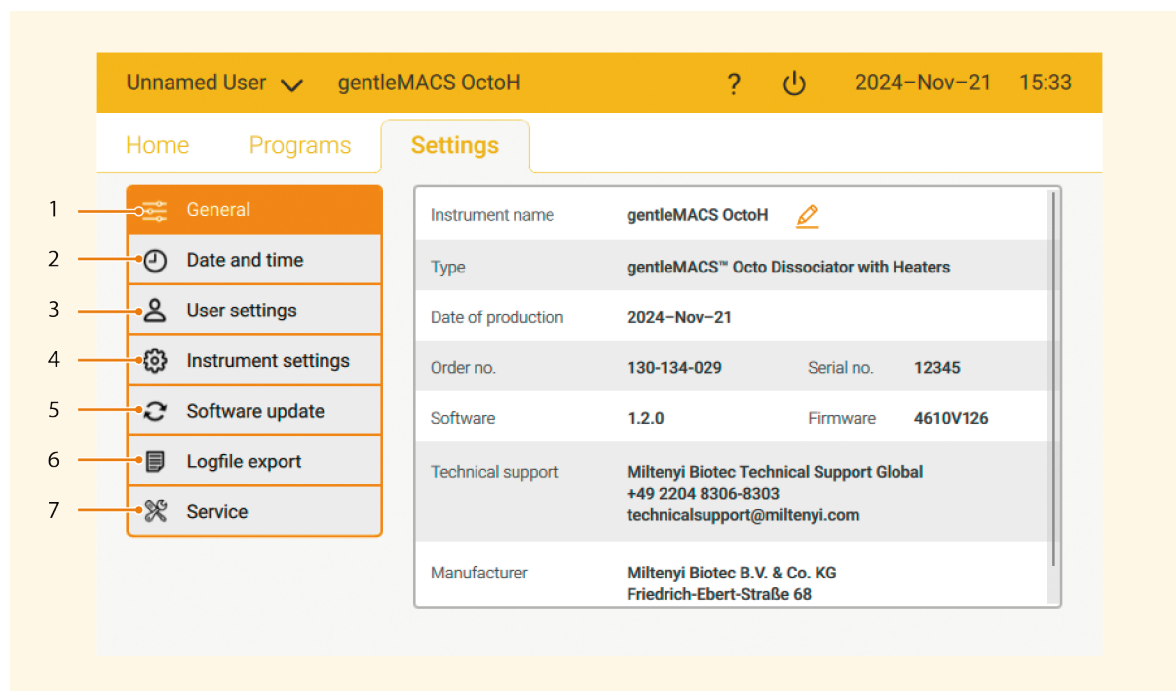
Figure 4.4: Overview Settings hub

- | | | |
|---|--------------------------------|--|
| 1 | General tab | General information about the instrument; tap the pen icon to change the instrument name |
| 2 | Date and time tab | Set date and time |
| 3 | User settings tab | Overview of all user profiles on this instrument, option to switch users, delete, clone, or edit user profiles, and to add new user profiles |
| 4 | Instrument settings tab | Settings for notifications, LED feedback, and display brightness |
| 5 | Software update tab | Check and update the software version |
| 6 | Logfile export tab | Export log files to a USB flash drive, see Exporting log files on page 78 |
| 7 | Service tab | Access to service mode for authorized service personnel |

4.6.1 Editing instrument name

- 1 Go to **Settings > General**.
- 2 Tap on the pen symbol next to the instrument name.
- 3 Edit the instrument name.
- 4 Confirm the changes.

4.6.2 Editing date and time

- 1 Go to **Settings > Date and time**.
- 2 Tap on the number to be changed. The selected numbers turn orange.
- 3 Edit date and time.
- 4 To confirm the changes, tap **Apply**.

4.6.3 User settings

Settings for notifications and brightness, favorite programs and the date of last use of programs by a user are saved in the profile of the signed-in user. User profiles help the different users to access their frequently used programs easier.

Adding new user profiles

- 1 Go to **Settings > User settings**.
- 2 Tap **New**.
- 3 Enter the user profile name.
- 4 To confirm, tap enter.
The new user profile is now included in the user profile list.

Switching user profiles

To switch between user profiles, do one of the following:

- Tap on the user name in the upper left corner of the top indicator bar and select the user profile to use.
- Go to **Settings > User settings** and select the user profile to use. Tap **Switch** to change the user profiles. The person icon indicates the currently active user.

Editing user profile names

- 1 Go to **Settings > User settings**.
- 2 Tap on the user profile to be edited.
The selected user profile is shown in orange.
- 3 Tap **Edit**.
- 4 Change the user profile name.
- 5 To confirm, tap enter.

Copying user profiles

Copying user profiles creates a copy of an existing user profile with the same settings defined in the user profile.

- 1 Go to **Settings > User settings**.
- 2 Tap on the user profile to be copied.
The selected user profile is shown in orange.
- 3 Tap **Clone**.
- 4 Enter a name for the new user profile.
- 5 To confirm, tap enter.

Deleting user profiles

The currently active user profile cannot be deleted. The user profile **Unnamed User** is the default user and cannot be deleted or edited.

- 1 Go to **Settings > User settings**.
- 2 Tap on the user profile to be deleted.
The selected user profile is shown in orange.
- 3 Tap **Delete**.
A pop-up window appears.
- 4 To confirm deletion of the user profile, tap **Delete**.

4.6.4 Instrument settings

Notifications and LED feedback for errors cannot be disabled.

The user can change the following settings for the respective user profile:

- information notifications,
- break state notifications,
- LED feedback of the instrument, and
- display brightness.

Tap the respective toggle switch to enable or disable the notifications or LED feedback. To receive further information about the respective setting, tap on the info icon.

Move the slider to increase or decrease the display brightness.

4.6.5 Software update

New software versions can be installed from a USB flash drive by the user. Make sure that the USB flash drive is formatted as exFAT for compatibility with the operating system. If required, change the format on a PC.

To update the Miltenyi Biotec program database and import new Miltenyi Biotec programs, see **Importing programs on page 65**.

- 1 Insert the USB flash drive to the USB port of the gentleMACS Octo Dissociator with Heaters.
- 2 Go to **Settings > Software update**.
- 3 Tap **Install**.
- 4 Choose the new software version to be installed from the USB flash drive.
- 5 Confirm to resume installation.
The instrument system checks if the software version is compatible with the instrument and installs the software.
- 6 The instrument reboots automatically with the newly installed software version and notifies the user about successful installation.

4.7 LEDs

Various statuses of the instrument are indicated by LEDs illuminating the Safety Shield and the gentleMACS Tubes. The LED feedback can be turned off, except for error feedback. See **User settings on page 67**.

LED status	Instrument status
Blue, solid	all positions in status running
Green, solid	at least one position is in status free , partly or ready
Green, blinking	at least one position is in status done
Blue, blinking	at least one position in status paused or break
Red, blinking	an error occurred. See Troubleshooting table on page 77 .

Table 4.3: Color code of the status LEDs

5

Performing runs

5.1 Running a single sample

- 1 Attach a tube upside down to the Sleeve of the respective position on the instrument. The instrument automatically detects the inserted tube and selects the position.

Insert and remove tubes in an upright position without tilting. Tilting might result in damage of the tube seal and subsequent leakage. Ensure that the tube is securely fastened to its Sleeve and socket. It must 'click' into position. Do not force the tube into the position. In case of resistance slightly turn the tube clockwise to find the right position to click in.

- 2 (Optional) If running a program including an incubation step, apply a gentleMACS Heater to the respective tube position. Slightly open the clamp of the gentleMACS Heater by pressing the upper handles and adjust it over the tube. Make sure the lower side of the gentleMACS Heater is facing towards you therefore allowing the electrical contact to close tightly. The instrument automatically detects the attached gentleMACS Heater.

Make sure that the gentleMACS Heater is securely fastened around the tube and to the socket. Do not use gentleMACS Heaters without a tube.

- 3 Tap on the desired program to run in the programs list.
The selected program is applied to the selected position box. The position status changes from **partly** to **ready**.
- 4 Tap the local start symbol in the position box to start the run.
The progress bar and timer indicate the remaining time until the program is done.
- 5 On completion of the program, the status of the position box changes to **done**.
- 6 Remove the tube from the tube position.
The status of the tube position changes to **free**.

5.2 Running multiple samples simultaneously

Up to eight samples can be run simultaneously using the same or different programs.

In this example, eight samples will be processed using the same program without a heating step.

- 1 Tap **Select all** to select all free position boxes in the **Home** hub.
- 2 Tap on the desired program to run in the programs list. The program is applied to all selected position boxes.
- 3 Attach tubes upside down to the Sleeves of the respective positions on the instrument. The instrument automatically detects the inserted tubes.

Insert and remove tubes in an upright position without tilting. Tilting might result in damage of the tube seal and subsequent leakage. Ensure that the tube is securely fastened to its Sleeve and socket. It must 'click' into position. Do not force the tube into the position. In case of resistance slightly turn the tube clockwise to find the right position to click in.

- 4 Press the **Start** button to start the run on all positions in the status **ready**.
The progress bar and timer indicate the time remaining for each position box.
- 5 On completion of the program, the status of the position boxes changes to **done**.

For further processing, use the **New** button to reset the position boxes from the **done** status into the **partly** status to assign a new program.

- 6 Remove the tubes from the tube positions.
The status of the position boxes changes to **free**.

5.3 Running a perfusion program

Before running a perfusion program, install the gentleMACS Perfusion Sleeves. For more information on the gentleMACS Perfusers and gentleMACS Perfusion Sleeves, refer to the respective data sheet.

- 1 Remove the gentleMACS Sleeve from the position you want to run a perfusion program.
To remove the Sleeve, pull it straight up.
- 2 Place the gentleMACS Perfusion Sleeve on the vacant metal pin.
- 3 Install the Sleeve by pressing it down completely.
- 4 After installation, the spring must be relaxed again.
- 5 Repeat step 1–4 for each position you want to run a perfusion program on.
- 6 Obey the steps described in the respective data sheet of the respective MACS Perfusion kit used.

For perfusion programs, manual steps must be performed during the **pause** status of the run.

5.4 Pausing and resuming samples

To pause a running sample, do one of the following:

- Tap the local **Pause** button of the respective position box. The status of the position changes from **running** to **paused**.
- Tap the global **Pause** button. All running samples will be paused. The status of all affected positions changes from **running** to **paused**.

To resume a paused sample, do one of the following:

- Tap the local **Start** button of the respective position box. The status of the position changes from **paused** to **running**.
- Tap the global **Start** button. All running samples will be paused. The status of all affected positions changes from **paused** to **running**.

To resume a break, tap the local **Start** button of the respective position box.

5.5 Aborting a running sample

- 1 Tap on the position boxes with runs to abort.
- 2 Tap **Abort**.
- 3 Confirm the notification by tapping **Abort**. The status of the selected position boxes changes from **running** to **done**.

6

Maintenance

WARNING

Contamination or infection could result in death or serious injury depending on the material used.

- Wear personal protective equipment (such as gloves, safety glasses, etc.) as indicated in the safety data sheet for the particular substance.

WARNING

Electric shock, short circuit, overheating, fire, and explosion could result in death or serious injury.

- Unplug the instrument before cleaning it.
- Use only small amounts of cleaning agents on a soft cloth to wipe the instrument. Do not spray or pour liquid cleaning agents onto or into the instrument.

Larger amounts of accidentally spilled liquid are automatically collected in the Trough of the gentleMACS Octo Dissociator with Heaters. Use a pipette to aspirate the liquid before removing the Trough for decontamination and cleaning.

Instrument parts wear over time, especially Sleeves. Depending on the sample material and frequency of use, these parts require regular maintenance and/or replacement. Contact Miltenyi Biotec Technical Support regarding spare parts. Miltenyi Biotec also offers service contracts for the gentleMACS Octo Dissociator with Heaters. Visit www.miltenyibiotec.com/gentlemacs-service for further information.

6.1 Decontamination and cleaning

In case of spillage or contamination, clean the instrument appropriately.

For disinfection, wipe the instrument with 70% ethanol or 70% isopropanol.

Clean the instrument only with cleaning agents that are noncorrosive and not damaging the plastic housing. Suitable cleaning agents include household detergents, 70% ethanol, and 70% isopropanol. Do not use acetone or any aggressive agents. If in doubt, if a certain cleaning agent is suitable for disinfection and/or cleaning the gentleMACS Octo Dissociator with Heaters, contact Miltenyi Biotec Technical Support.

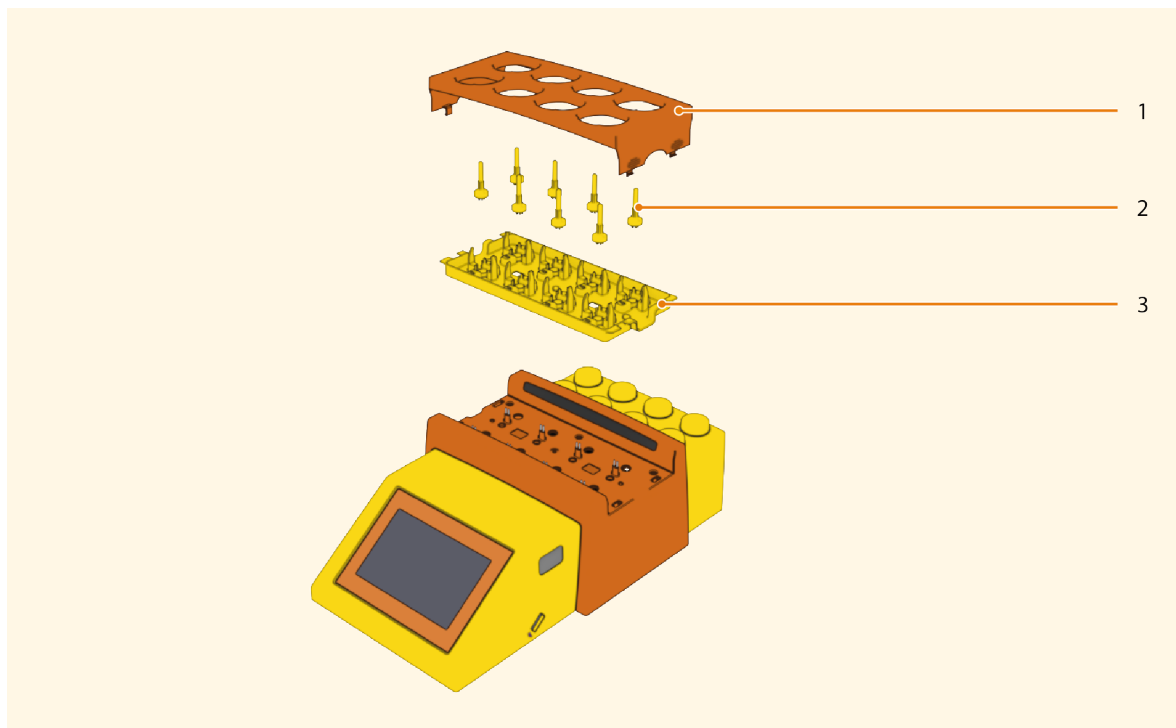
Figure 6.1: Steps 2-4 of the gentleMACS Octo Dissociator with Heaters disassembly for cleaning

- 1** Unplug the power cable from the electrical outlet before cleaning the instrument.
- 2** Remove the Safety Shield (1).
- 3** Remove the Sleeves from the Sleeve holders (2).
- 4** Remove the Trough from the baseplate (3).
To remove the Trough, pull on one side. Do not pull on both sides at the same time.
- 5** Carefully wipe the surface of the instrument and all removed parts with a cloth soaked with a detergent-containing aqueous solution, 70% ethanol or 70% isopropanol.
- 6** Repeat the procedure with another cloth soaked with distilled water.
- 7** Reassemble the removed parts. Start with the Trough, followed by the Sleeves, and the Safety Shield. Make sure that the orientation of the Trough and Safety Shield is correct.

7

Troubleshooting

7.1 Troubleshooting table

Issue	Possible reason	Solution
Liquid droplets appear in the Trough.	Tube is not closed properly.	Close the tube tightly beyond first resistance.
	Sample is jammed between the cap and the tube and damages the seal. This might cause liquid to run through the seal.	The top area of the cap and the tube must be perfectly clean for proper functioning of the seal.
	Tube seal is leaky.	Replace the tube.
	Sleeve is worn.	Exchange Sleeve.
Tube is blocked during a run.	Inappropriate sample material.	Check the respective data sheet for technical specifications of the tubes.
Program is automatically aborted and the instrument displays the message "Motor is blocked. Program was aborted."	If the motor is blocked due to a high resistance of the sample material, the instrument will try three times to get started again	Process a smaller amount of tissue, or use a different gentleMACS Program.
	If the blockage can be resolved, the instrument will continue the run of the program. If the blockage cannot be resolved, the program will automatically be aborted.	Check if tissue is trapped between the stator and rotor. Release tissue, reinstall tube, and tap Continue to resume the program.
Tube comes off the instrument.	Tube is not correctly mounted into the snap-fit of the Trough.	Ensure correct installation of the tube.
	Tube is not closed properly.	Close the tube tightly beyond first resistance.
A warning message appears in a pop-up window when saving an edited program. "Input error - Maximum rounds per run (3200) exceeded!"	The duration of the program exceeds 3,200 rounds per run (rpr). It is not recommended to run programs exceeding this limitation! Miltenyi Biotec does not guarantee proper function of the instruments and disposables above this value.	Press OK to save the program. Press Abort to re-edit the program.

Issue	Possible reason	Solution
New program containing iterations cannot be saved. The instrument displays an error message with the cause of the error.	"Loop command is faulty. End loop command is missing." indicates that the end loop command has not been specified.	Close iterations by adding an end loop command.
	"Loop command is faulty. The command is nested." indicates that there are more than one loop commands.	Use only one loop command and close the iteration by adding an end loop command
	"Loop command is faulty. Loop has no body." indicates that there is no command between the start and end loop command.	At least one program step must be included between a start and end loop command. Add steps and define commands that shall be iterated.
A warning message "The heater was removed!" appears in a pop-up window while running a heaters program.	The heater was removed.	Reinstall heater.
	The heater is not installed properly.	Make sure that the heater is installed properly. Carefully clean the heater contacts of the instruments and the resp. heater using a cloak soaked with water or 70% ethanol.
	The electrical contacts of the instrument or the heater are dirty or corroded.	Carefully check the heater for signs of damage. If no visible signs of damage can be detected, switch the instrument off and on again and retry to run the program.
		Contact Miltenyi Biotec Technical Support.
Software for updating the Miltenyi Biotec program database is not detected.	The Software update tab is used for updating the Miltenyi Biotec program database.	Updating the Miltenyi Biotec program database is not a software update. To import new Miltenyi Biotec programs, go to the Programs hub and tap Import. See Importing programs on page 65.
USB flash drive is not recognized.	The USB flash drive has the wrong format.	Format the USB flash drive as exFAT.

Table 7.1: Troubleshooting for the gentleMACS Octo Dissociator with Heaters

7.2 Exporting log files

Log files can help Miltenyi Biotec Technical Support with troubleshooting. Log files can be exported to a USB flash drive. Make sure that the USB flash drive is formatted as exFAT for compatibility with the operating system. If required, change the format on a PC.

- 1 Insert the USB flash drive to the USB port of the gentleMACS Octo Dissociator with Heaters.
- 2 Go to **Settings > Logfile export**
- 3 Tab **Export logfile** to start the export.
- 4 After notification that the export is finished, remove the USB flash drive from the instrument.

8

Technical support

For technical support, contact your local Miltenyi Biotec representative or Miltenyi Biotec Technical Support at Miltenyi Biotec headquarters:

Miltenyi Biotec B.V. & Co. KG
Friedrich-Ebert-Straße 68
51429 Bergisch Gladbach
Germany
Phone: +49 2204 8306 3803
technicalsupport@miltenyi.com

Visit **www.miltenyibiotec.com** for local Miltenyi Biotec Technical Support contact information.

Technical data and specifications

Technical data	Specification
Colors	Yellow / orange
Dimensions (width x depth)	285 mm × 496 mm
Height	188.5 mm (without heaters) 245 mm (with heaters)
Weight	11.8 kg
Main power	
Power consumption	450 W
Supply voltage	100–240 V~, 50/60 Hz
Fuses	2 x 3.15 AH, 250 V
Motor	
Rotational speed	20–4,000 rpm
Rotational direction	Clockwise or counterclockwise
Rotational acceleration	100 r/s ²
Torque	Max. 150 mNm; Overload restriction over the range of 200–4,000 rpm
Conditions of operation	
Working temperature without use of heaters	+5 °C to +35 °C
Working temperature with use of heaters	+18 °C to +30 °C
Humidity	15–85%, relative, noncondensing
Altitude	Max. 2,000 m
Storage temperature	-15 °C to +70 °C
Ports	1 × USB 2.0 1 × Ethernet (10/100/1000 MBit/s) 1 × CAN (1 MBit/s)
RAM	4 GB DDR3L
Mass storage	32 GB mSATA SSD
Mainboard battery	Button cell, lithium metal
Monitor	7" TFT LCD touchscreen
Maximum number of programs	500
Number of heatable positions	8

Technical data	Specification
Maximum noise emission	
Sound pressure level at workstation	85 dB(A) +/- 2.5 dB(A)
Sound power level	91 dB
Measurement operation conditions: 8x C Tubes containing 500 µL of liquid, processed at 4,000 rpm	
Working temperature heaters	37 °C

Table 9.1: Technical data of the instrument

The gentleMACS™ Octo Dissociator with Heaters is classified as protection class I device and must be plugged into a grounded electrical outlet.

Note that for USA and Canada, the main power supply cord and plug of the instrument shall comply with the following specifications: UL listed and KAM cord, minimum type SJ, minimum 18 AWG, 3 conductors. Rated for a minimum temperature of 60 °C. Provided with grounding-type (NEMA 5-15P) attachment plug, rated 125 VAC, 10 A. Opposite end terminates in IEC 320-style connector, rated 125 VAC, 10A.

Supply voltage fluctuations up to +/- 10% of the nominal voltage.

Transient overvoltages present on the mains supply: category II.

The instrument is suitable for pollution degree 2 environments.

The instrument has been designed for a lifetime of 10 years.

9.1 FCC compliance statement

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

Changes or modifications made to this equipment not expressly approved by the party responsible for compliance may void the FCC authorization to operate this equipment. For other safety considerations, refer to the product label or visit www.miltenyibiotec.com. Design and specifications are subject to change without notice.

9.2 EC/EU Declaration of Conformity

This declaration of conformity is issued under the sole responsibility of the manufacturer:

Miltenyi Biotec B.V. & Co. KG
Friedrich-Ebert-Straße 68
51429 Bergisch Gladbach
Germany

This declaration relates exclusively to the machinery in the state in which it was placed on the market, and excludes components which are added and/or operations carried out subsequently by the final user.

The declaration of conformity refers to the machinery identified as follows:

Description: Laboratory equipment
Model: gentleMACS Octo Dissociator with Heaters (A44109)

The machinery complies with all essential requirements of the following directives:

2006/42/EC Machinery
2011/65/EU Restriction of the use of certain hazardous substances in electrical & electronic equipment
2014/30/EU Electromagnetic compatibility

The machinery is in conformity with the following harmonized standards:

EN 61010-1:2010
EN 61010-2-010:2014
EN 61010-2-051:2015
EN 61326-1:2013

Person authorized to compile the relevant technical documentation:

Dr. Bernd Schröder
Global Head Regulatory Affairs
Miltenyi Biotec B.V. & Co. KG
Friedrich-Ebert-Straße 68
51429 Bergisch Gladbach
Germany

9.3 UK Declaration of Conformity

This declaration of conformity is issued under the sole responsibility of the manufacturer:

Miltenyi Biotec B.V. & Co. KG
Friedrich-Ebert-Straße 68
51429 Bergisch Gladbach
Germany

This declaration relates exclusively to the machinery in the state in which it was placed on the market, and excludes components which are added and/or operations carried out subsequently by the final user.

The declaration of conformity refers to the machinery identified as follows:

Description: Laboratory equipment
Model: gentleMACS Octo Dissociator with Heaters (A44109)

The machinery complies with all essential requirements of the following legislations:

Supply of machinery (Safety) regulations 2008
The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012
Electromagnetic Compatibility Regulations 2016

The machinery is in conformity with the following UK designated standards:

EN 61010-1:2010
EN 61010-2-010:2014
EN 61010-2-051:2015
EN 61326-1:2013

Person authorized to compile the relevant technical documentation:

Dr. Bernd Schröder
Global Head Regulatory Affairs
Miltenyi Biotec B.V. & Co. KG
Friedrich-Ebert-Straße 68
51429 Bergisch Gladbach
Germany

10

Legal notes

10.1 Limited warranty

Except as stated in a specific warranty statement which may accompany this product or as otherwise agreed in writing by an authorized representative of Miltenyi Biotec, Miltenyi Biotec's warranty to you, the original purchaser and end user ("you" or "your"), with respect to the product accompanied by this limited warranty shall be subject to the following provisions and the general terms and conditions of sale of the company within the Miltenyi Biotec group which supplied the product in effect at the date of purchase. Those terms and conditions of sale may vary by country and region. Nothing in this document should be construed as constituting an additional warranty.

Miltenyi Biotec warrants that this product will operate or perform substantially in conformance with Miltenyi Biotec's published specifications and be free from material defects in material and workmanship, when subjected to normal, proper, and intended usage by properly trained personnel, for the period of time set forth in the product documentation or package inserts accompanying the product (the "Warranty Period").

Miltenyi Biotec agrees, during the Warranty Period, to repair or replace, at Miltenyi Biotec's option, the defective product so as to cause the same to operate in substantial conformance with said published specifications; provided that you shall (a) promptly notify Miltenyi Biotec in writing upon the discovery of any nonconformity or defect, which notice shall include the product model and serial number (if applicable) and details of the warranty claim; and (b) return the nonconforming or defective product to Miltenyi Biotec, freight prepaid, only after receipt of a Return Material Authorization ("RMA") from Miltenyi Biotec, which may include biohazard decontamination procedures and other product-specific handling instructions, if applicable.

Miltenyi Biotec shall have no obligation to make repairs, replacements, or corrections to the product or any component thereof required, in whole or in part, as the result of (i) normal wear and tear, (ii) improper handling, installation, operation, storage, service, maintenance, or repair, (iii) failure to follow the instructions, cautions, warnings, and notes set forth in the product documentation provided with the product or provided by Miltenyi Biotec from time to time, (iv) abnormal use, misuse, neglect, abuse, mishandling, misapplication, modification, or alteration of the product, (v) use of the product in a manner for which it was not designed, (vi) causes external to the product such as, but not limited to, power failure or electrical power surges, (vii) use of the product in combination with equipment, accessories, consumables, or software not supplied or approved by Miltenyi Biotec, or (viii) accident, disaster, or acts of God. ANY INSTALLATION, MAINTENANCE, REPAIR, SERVICE, OR ALTERATION TO OR OF, OR OTHER TAMPERING WITH, THE PRODUCT PERFORMED BY ANY PERSON OR ENTITY OTHER THAN MILTENYI BIOTEC AUTHORIZED PERSONNEL WITHOUT MILTENYI BIOTEC'S PRIOR WRITTEN APPROVAL, OR ANY USE OF REPLACEMENT PARTS NOT SUPPLIED BY MILTENYI BIOTEC, SHALL IMMEDIATELY VOID AND CANCEL ALL WARRANTIES WITH RESPECT TO THE AFFECTED PRODUCT.

Miltenyi Biotec's warranty does not cover products sold AS IS or WITH ALL FAULTS, or which had its serial number defaced, altered, or removed, or any consumables or parts identified as being supplied by a third party.

Miltenyi Biotec must be informed promptly if a claim is made under this limited warranty. If a material or manufacturing defect occurs within the Warranty Period, Miltenyi Biotec will take the appropriate steps, at Miltenyi Biotec's option, to make repairs, replacements, or corrections to the product or any component thereof as may be required to restore the full usability of your product. If Miltenyi Biotec determines that a product for which you have requested warranty services is not covered by the warranty hereunder, you shall pay or reimburse Miltenyi Biotec for all costs of investigating and responding to such request at Miltenyi Biotec's then prevailing time and materials rates. If Miltenyi Biotec provides repair services or replacement parts that are not covered by this warranty, you shall pay Miltenyi Biotec therefor at Miltenyi Biotec's then prevailing time and materials rates.

THE OBLIGATIONS CREATED BY THIS WARRANTY STATEMENT TO REPAIR OR REPLACE A DEFECTIVE PRODUCT ARE EXCLUSIVE AND SHALL BE THE SOLE REMEDY OF BUYER IN THE EVENT OF A DEFECTIVE PRODUCT. EXCEPT AS EXPRESSLY PROVIDED IN THIS WARRANTY STATEMENT, MILTENYI BIOTEC HEREBY DISCLAIMS ALL OTHER WARRANTIES, WHETHER EXPRESS OR IMPLIED, ORAL OR WRITTEN, WITH RESPECT TO THE PRODUCT, INCLUDING WITHOUT LIMITATION ALL IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS FOR ANY PARTICULAR PURPOSE. MILTENYI BIOTEC DOES NOT WARRANT THAT THE PRODUCT IS ERROR-FREE OR WILL ACCOMPLISH ANY PARTICULAR RESULT.

UNDER NO CIRCUMSTANCES WILL MILTENYI BIOTEC BE LIABLE FOR ANY LOSS OF USE, INTERRUPTION OF BUSINESS, OR ANY INDIRECT, SPECIAL, INCIDENTAL, PUNITIVE, OR CONSEQUENTIAL DAMAGES OF ANY KIND (INCLUDING LOST PROFITS) REGARDLESS OF THE FORM OF ACTION WHETHER IN CONTRACT, TORT (INCLUDING NEGLIGENCE), STRICT PRODUCT LIABILITY, OR OTHERWISE, EVEN IF MILTENYI BIOTEC HAS BEEN ADVISED OF THE POSSIBILITY OF SUCH DAMAGES. IN NO EVENT SHALL THE TOTAL LIABILITY OF MILTENYI BIOTEC HEREUNDER EXCEED THE GREATER OF 100.00 EUROS OR THE AMOUNT YOU ACTUALLY PAID FOR THE PRODUCT GIVING RISE TO SUCH LIABILITY, REGARDLESS OF THE CAUSE OF ACTION, IN CONTRACT, TORT, STRICT LIABILITY, OR OTHERWISE.

NOT ALL JURISDICTIONS ALLOW SUCH LIMITATIONS OF DAMAGES SO THE FOREGOING LIMITATIONS MAY NOT APPLY TO YOU. This warranty statement gives you specific legal rights and you may have other rights, which may vary from jurisdiction to jurisdiction.

10.2 Trademarks

gentleMACS, MACS, and the Miltenyi Biotec logo are registered trademarks or trademarks of Miltenyi Biotec B.V. & Co. KG and/or its affiliates in various countries worldwide.

11

Index

D

Disassembly **76**

E

Editing

Date and time **66-67**

Program **64**

User profile **67**

G

gentleMACS Heaters **52**

gentleMACS Octo Coolers **53**

gentleMACS Tubes **51**

H

Help **60**

L

Logfile export **66, 78**

P

Program

Create **63**

Delete **65**

Edit **64**

Export **65**

Filter **64**

Import **65**

S

Safety Shield **49, 58**

Sample

Abort **73**

Pause **72**

Run **71**

Sleeve **49, 58**

Status

LED **69**

Position boxes **62**

Switching off **59**

Switching on **59**

T

Trough **49, 58**

U

Update

Miltenyi Biotec program database **65**

Software **66, 68**

User profile

Add **67**

Copy **67**

Delete **68**

Edit **67**

Switch **67**

Miltenyi Biotec

Germany/Austria

Miltenyi Biotec B.V. & Co. KG
Friedrich-Ebert-Straße 68
51429 Bergisch Gladbach
Germany
Phone +49 2204 8306-0
Fax +49 2204 85197
macsde@miltenyi.com

USA/Canada

Miltenyi Biotec, Inc.
1201 Clopper Road
Gaithersburg, MD 20878
USA
Phone 800 FOR MACS
Phone +1 866 811 4466
Fax +1 877 591 1060
macsus@miltenyi.com

Australia

Miltenyi Biotec Australia Pty. Ltd.
Unit 11, 2 Eden Park Drive
Macquarie Park NSW 2113
Australia
Phone +61 2 8877 7400
Fax +61 2 9889 5044
macsau@miltenyi.com

Benelux

Miltenyi Biotec B.V.
Dellaertweg 9C
2316 WZ Leiden
The Netherlands
macsnl@miltenyi.com
**Customer service for:
The Netherlands**
Phone 0800 4020120
Fax 0800 4020100

Belgium

Phone 0800 94016
Fax 0800 99626

Luxembourg

Phone 800 24971
Fax 800 24984

China

Miltenyi Biotec Technology &
Trading (Shanghai) Co., Ltd.
Room A401, 4/F
No. 1077, Zhangheng Road
Pudong New Area
201203 Shanghai, P.R. China
Phone +86 21 6084 0210
Fax +86 21 6235 0953
macscn@miltenyi.com.cn

France

Miltenyi Biotec SAS
10 rue Mercoeur
75011 Paris, France
Phone +33 1 56 98 16 16
macsfr@miltenyi.com

Hong Kong

Miltenyi Biotec Hong Kong Ltd.
Unit 301, Lakeside 1
No. 8 Science Park West Avenue
Hong Kong Science Park
Pak Shek Kok, New Territories
Hong Kong
Phone +852 3751 6698
Fax +852 3619 5772
macshk@miltenyi.com.hk

India

Miltenyi Biotec India Pvt. Ltd.
Ground Floor, Lab No. E 101/1,
Building no. 3600 | Plot No. 4,
Sy. no. 101, Neovantage Park,
Synergy Square 3 Genome
Valley, Lalgadimalakpet
Medchal Malkajgiri, Shamirpet
(M) | Telangana 500101, India
Phone +91 040 45175910
macsin@miltenyi.com

Italy

Miltenyi Biotec S.r.l.
Via Paolo Nanni Costa, 30
40133 Bologna, Italy
Phone +39 051 6 460 411
Fax +39 051 6 460 499
macsit@miltenyi.com

Japan

Miltenyi Biotec K.K.
NEX-Eitai Building 5F
16-10 Fuyuki, Koto-ku,
Tokyo 135-0041, Japan
Phone +81 3 5646 8910
Fax +81 3 5646 8911
macsjp@miltenyi.com

Nordics and Baltics

Miltenyi Biotec Norden AB
Medicon Village
Scheeleorget 1
223 81 Lund, Sweden
macsse@miltenyi.com

Customer service for:

Sweden

Phone 0200 111 800
Fax +46 280 72 99

Denmark

Phone 80 20 30 10
Fax +46 46 280 72 99

Norway, Finland, Iceland, and Baltic countries

Phone +46 46 280 72 80
Fax +46 46 280 72 99

Singapore

Miltenyi Biotec Asia Pacific Pte Ltd.
438B Alexandra Road, Block B
Alexandra Technopark
#06-01
Singapore 119968
Phone +65 6238 8183
Fax +65 6238 0302
macssg@miltenyi.com

South Korea

Miltenyi Biotec Korea Co., Ltd.
Donggeuk 7F,
562 Nonhyeon-ro
Gangnam-gu
Seoul 06136, South Korea
Phone +82 2 555 1988
Fax +82 2 555 8890
macskr@miltenyi.com

Spain

Miltenyi Biotec S.L.
C/Virgilio 2, Edificio II, Planta -1
28223 Pozuelo de Alarcón, Madrid
Spain
Phone +34 91 512 12 90
Fax +34 91 512 12 91
macses@miltenyi.com

Switzerland

Miltenyi Biotec Swiss AG
Soodstrasse 52
8134 Adliswil
Switzerland
Phone +41 32 623 08 47
Fax +49 2204 85197
macsch@miltenyi.com

United Kingdom

Miltenyi Biotec Ltd.
Almac House, Church Lane
Bisley, Surrey GU24 9DR, UK
Phone +44 1483 799 800
Fax +44 1483 799 811
macsuk@miltenyi.com

www.miltenyibiotec.com

Miltenyi Biotec provides products and services worldwide. Visit www.miltenyibiotec.com/local to find your nearest Miltenyi Biotec contact.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research use only and not for therapeutic or diagnostic use.